

LAB DERATING INSTRUCTIONS SIMDOS® 02 DIAPHRAGM METERING PUMP



KNF Flodos AG Wassermatte 2 6210 Sursee, Switzerland Tel +41 (0)41 925 00 25 Fax +41 (0)41 925 00 35 www.knf.com

Contents	6	Page
1	General points	5
1.1	Information on the instructions	5
1.2	Warnings	6
1.3	Symbols	7
1.4	Disclaimer of liability	8
1.5	Manufacturer's address	8
1.6	Year of manufacture	8
1.7	Other applicable documents	8
1.8	Copyright	8
1.9	Warranty conditions	8
2	Safety	9
2.1	Intended use	9
2.2	Reasonably foreseeable misuse	9
2.3	Owner's responsibility	10
2.4	Personnel requirements	10
2.5	Product-specific dangers	11
2.6	Personal protective equipment	12
2.7	Safety equipment	13
2.8	Environmental protection considerations	13
2.9	Danger areas	13
2.10	Declaration of conformity	14
3	Technical data	15
3.1	Items included in delivery	15
3.2	Storage conditions	15
3.3	Pump head materials	15
3.4	Product key	16
3.5	Dimensions	17
3.6	Installation location requirements	18
3.7	Electrical connections and performance data	18
3.8	Remote control (RC version only)	19
3.9	Remote control via RS 232 (RC Plus version only)	19
	Flow rates	20
	Other parameters	20
	Hydraulic connections	20
	Transferred medium	21
3.14	Accuracy/ reproducibility	22
4	Assembly and function	
4.1	Metering pump structure	24
4.2	Operating principle	24
4.3	Valves	25
4.4	Pump diaphragm	27
4.5	Pump head	27
4.6	Operating controls	27
4.7	Operating modes	31
5	Shipment	
5.1	Checking delivery	32
5.2	Registering a complaint	32
5.3	Arranging for return	32
6	Installation and initial start-up	
6.1	Safety	33
6.2	Installation	34
6.3	System setup	37
6.4	Initial start-up	39
7	Operation	
7.1	Safety	40
7.2	Switching pump on / off	41
7.3	Entering settings	42
7.4	Setting the operating mode	43
7.5	Priming the pump	44

7.6	Starting the pump	44
7.7	Interrupting a transferring operation	44
7.8	Stopping the transferring operation	45
7.9	Transferring fluid	46
7.10	Setting metering operation	46
7.11	Setting the fluid type	47
7.12	To calibrate the pump	48
7.13	Cycle metering mode	49
7.14	Setting the system	49
7.15	Back to main menu	49
7.16	Select language setting	50
7.17 7.18	Activate autostart Set contrast	50 50
	Resetting the pump (Reset)	50
	Shutting down the pump in an emergency	51
	RC version (remote control)	
o 1 8.1	Remote control analog input	52 53
8.2	Digital input 1, Start / Stop	55
8.3	Digital input 2 Reset / Prime / Foot Switch	57
8.4	Digital output	59
	•	
9	RC Plus version (serial communication)	60
	RC Plus version (serial communication)	
10	Cleaning and maintenance	61
10 10.1	Cleaning and maintenance Safety	61 61
10 10.1 10.2	Cleaning and maintenance Safety Cleaning the pump	61 61 62
10 10.1 10.2 10.3	Cleaning and maintenance Safety Cleaning the pump Maintenance plan	61 61 62 63
10 10.1 10.2 10.3 10.4	Cleaning and maintenance Safety Cleaning the pump Maintenance plan Cleaning / replacing valve plates and pump diaphragm	61 61 62
10 10.1 10.2 10.3 10.4 10.5	Cleaning and maintenance Safety Cleaning the pump Maintenance plan Cleaning / replacing valve plates and pump diaphragm Checking that pump is leak-tight	61 61 62 63 64 70
10 10.1 10.2 10.3 10.4 10.5	Cleaning and maintenance Safety Cleaning the pump Maintenance plan Cleaning / replacing valve plates and pump diaphragm Checking that pump is leak-tight Shutdown	61 61 62 63 64 70
10 10.1 10.2 10.3 10.4 10.5 11 11.1	Cleaning and maintenance Safety Cleaning the pump Maintenance plan Cleaning / replacing valve plates and pump diaphragm Checking that pump is leak-tight Shutdown Safety	61 61 62 63 64 70 71 71
10 10.1 10.2 10.3 10.4 10.5 11 11.1	Cleaning and maintenance Safety Cleaning the pump Maintenance plan Cleaning / replacing valve plates and pump diaphragm Checking that pump is leak-tight Shutdown	61 61 62 63 64 70 71 71
10 10.1 10.2 10.3 10.4 10.5 11 11.1 12	Cleaning and maintenance Safety Cleaning the pump Maintenance plan Cleaning / replacing valve plates and pump diaphragm Checking that pump is leak-tight Shutdown Safety Troubleshooting Safety	61 61 62 63 64 70
10 10.1 10.2 10.3 10.4 10.5 11 12 12.1	Cleaning and maintenance Safety Cleaning the pump Maintenance plan Cleaning / replacing valve plates and pump diaphragm Checking that pump is leak-tight Shutdown Safety Froubleshooting	61 61 62 63 64 70
10 10.1 10.2 10.3 10.4 10.5 11 12 12.1 12.2 12.3	Cleaning and maintenance Safety Cleaning the pump Maintenance plan Cleaning / replacing valve plates and pump diaphragm Checking that pump is leak-tight Shutdown Safety Troubleshooting Safety Displaying of error messages Rectifying transfer problems	61 61 62 63 64 70 71 71 71 72 72 72 73
10 10.1 10.2 10.3 10.4 10.5 11 12 12.1 12.2 12.3	Cleaning and maintenance Safety Cleaning the pump Maintenance plan Cleaning / replacing valve plates and pump diaphragm Checking that pump is leak-tight Shutdown Safety Troubleshooting Safety Displaying of error messages	61 61 62 63 64 70 71 71 71 72 72 72 73
10 10.1 10.2 10.3 10.4 10.5 11 12.1 12.1 12.2 12.3 13 13.1	Cleaning and maintenance Safety Cleaning the pump Maintenance plan Cleaning / replacing valve plates and pump diaphragm Checking that pump is leak-tight Shutdown Safety Troubleshooting Safety Displaying of error messages Rectifying transfer problems Spare parts and accessories.	61 61 62 63 64 70 71 71 71 72 72 72 73 74
10 10.1 10.2 10.3 10.4 10.5 11 12 12.1 12.2 12.3 13 13 13.2 13.3	Cleaning and maintenance Safety Cleaning the pump Maintenance plan Cleaning / replacing valve plates and pump diaphragm Checking that pump is leak-tight Shutdown Safety Troubleshooting Safety Displaying of error messages Rectifying transfer problems Spare parts and accessories	61 61 62 63 64 70 71 71 71 72 72 72 73 73 74 74
10 10.1 10.2 10.3 10.4 10.5 11 12 12.1 12.2 12.3 13 13 13.1 13.2 13.3 13.4	Cleaning and maintenance Safety Cleaning the pump Maintenance plan Cleaning / replacing valve plates and pump diaphragm Checking that pump is leak-tight Shutdown Safety Troubleshooting Safety Displaying of error messages Rectifying transfer problems Spare parts and accessories Spare parts Hoses Pressure control valve Filter	61 61 62 63 64 70 71 71 71 72 72 72 73 73 74 74 74
10 10.1 10.2 10.3 10.4 10.5 11 12 12.1 12.2 12.3 13 13 13.1 13.2 13.3 13.4 13.5	Cleaning and maintenance Safety Cleaning the pump Maintenance plan Cleaning / replacing valve plates and pump diaphragm Checking that pump is leak-tight Shutdown Safety Troubleshooting Safety Displaying of error messages Rectifying transfer problems Spare parts and accessories Spare parts Hoses Pressure control valve Filter Foot switch	61 61 62 63 64 70 71 71 71 72 72 73 73 73 74 74 74 74 74 74 74
10 10.1 10.2 10.3 10.4 10.5 11 12 12.1 12.2 12.3 13 13.1 13.2 13.3 13.4 13.5 13.6	Cleaning and maintenance Safety Cleaning the pump Maintenance plan Cleaning / replacing valve plates and pump diaphragm Checking that pump is leak-tight Shutdown Safety Troubleshooting Safety Displaying of error messages Rectifying transfer problems Spare parts and accessories Spare parts Hoses Pressure control valve Filter	61 61 62 63 64 70 71 71 71 72 72 73 73 73 74 74 74 74 74 74

1 General points

1.1 Information on the instructions

Contents	The operating instructions contain important notes on how to use the pump. In order to ensure safe working and proper functioning it is essential to observe all the specified safety precautions.
Storage location	These operating instructions are part of the product, and must be stored in its immediate vicinity in a location accessible to personnel at all times.
Passing on	These operating instructions are part of the product, and must be passed on to the next owner if the device is resold.
Project pumps	Customer-specific project pumps (pump models which begin with "PL" or "PML") may differ from the operating instructions, in which case the agreed specification is also applicable. It is listed with the other applicable documents.
Illustrations in the instructions	Illustrations in these operating instructions may be to scale, but do not have be. The illustrations may differ slightly from the actual product.

1.2 Warnings

Warnings in the operating instructions are identified with the danger symbol, keywords and colours. These provide an indication of the extent of the danger.

A DANGER

Indicates a dangerous situation which may lead directly to death or serious injury if it is not avoided.

Indicates a dangerous situation which may lead to death or serious injury if it is not avoided.

A CAUTION

Indicates a dangerous situation which may lead to moderate or minor injuries if it is not avoided.

NOTE

Indicates a situation which may cause damage to property if it is not avoided.







1.3 Symbols

The following symbols appear in the operating instructions, on labels on the device, and on its packaging:



A WARNING

Danger of injuries and property damage caused by illegible labels

The labelling on the device may become illegible with time.

- \triangleright Maintain labels on the device in a legible condition.
- Replace any illegible labels. \geq

The following symbols in the operating instructions and on the device and its packaging indicate environmental protection considerations:



Recycling

Not to be disposed of in household waste

The following symbols in safety precautions and on the device indicate the nature of the danger:



General danger symbol Further details of the nature of the danger are provided in the warning highlighted by this symbol.



Electrical danger



Fire danger





Danger of automatic start-up

Danger of caustic burns or burns

This symbol calls for disconnection from the mains.

1.4 Disclaimer of liability

In preparing the contents of these operating instructions, account has been taken of applicable regulations.

The manufacturer can accept no liability for any damage or malfunctions caused by failure to follow the operating instructions.

The manufacturer can accept no liability for any damage or malfunctions caused by modification or conversion of the device or improper handling.

The manufacturer can accept no liability for any damage or malfunctions caused by the use of non-approved spare parts and accessories.

1.5 Manufacturer's address

KNF Flodos AG Wassermatte 2 6210 Sursee, Switzerland Tel +41 (0)41 925 00 25 Fax+41 (0)41 925 00 35 www.knf.com

1.6 Year of manufacture

The year of manufacture is shown on the device's type plate.

1.7 Other applicable documents

The documents listed must also be taken into consideration.

- Chemical resistance list
- Sales documents
- General Trade Conditions

1.8 Copyright

The information, texts and illustrations in these operating instructions are protected by copyright. The contents of these operating instructions must not be copied, translated or passed to third parties without the written consent of the manufacturer.

1.9 Warranty conditions

The applicable warranty conditions are set out in the General Terms and Conditions of Business and the sales documents.

The warranty does not cover the following:

- Malfunctions caused by particles/fibres in the valve system
- Sticking of the valve system as a result of inadequate flushing

2 Safety

2.1 Intended use

This pump is intended exclusively for use as follows:

- Transferring and metering liquids.
- Temporary operation in laboratories for research purposes.
- Operation in accordance with the operating parameters specified in the technical data (see Chapter 3) and other applicable documents.
- Transferring media which meet the requirements in the technical data and other applicable documents.
- Operation when fully assembled.
- Upright operation on a firm table top.

2.2 Reasonably foreseeable misuse

The pump must not be

- operated in an explosive atmosphere.
- used to transfer explosive media.
- used to produce foodstuffs or pharmacological products. Additional certifications are required for this purpose.
- used to transfer media whose compatibility with the pump head, valves and seals has not been demonstrated.

2.3 Owner's responsibility

The operator is responsible for compliance with the safety precautions in the present operating instructions and with any safety, accident prevention and environmental protection regulations which apply to the medium to be transferred and the range of applications of the device.

These include the following in particular:

- The operator must perform a risk assessment to identify any additional risks which may arise from the specific working conditions at the pump's installation location and the nature of the medium to be transferred. He/she must convert these into instructions for operating the device.
- The operator must ensure that any employees working with the pump have read the operating instructions.
- The operator must train his personnel at regular intervals and inform them of the risks involved.
- The operator must provide his personnel with the necessary protective equipment as specified in the safety data sheet for the medium to be transferred.
- The operator must maintain the pump in perfect working order.
- The operator must ensure compliance with the maintenance intervals described in the operating instructions.

2.4 Personnel requirements

A WARNING Danger of injuries and property damage caused by insufficiently qualified staff Improper use of the pump may result in serious injury and damage to property. Any activities must only be performed by adequately qualified personnel The following qualifications are specified for various activities in the operating instructions: Trained personnel These have been instructed in the task assigned to them and made aware of potential dangers and risks. This training is the minimum qualification for working on the device. Where no other qualification is specified, the activity must be performed by a trained individual as a minimum. The specialist training, knowledge and experience of such person-Specialist personnel nel, together with their knowledge of the relevant regulations, standards and directives, allow them to perform their assigned duties correctly, and to independently identify and avoid potential dangers and risks. Unauthorised personnel Individuals whose reactions have been impaired by the consumption of medicines, alcohol, drugs or by other causes and conditions are not authorised to perform activities on the device.

10

2.5 **Product-specific dangers**

This Chapter describes residual risks that have been identified in a risk assessment. Safety precautions and warnings in this Chapter and in other Chapters in the operating instructions must be observed in order to avoid dangerous situations.

Danger of injuries and property damage caused by leaks in connections and the pump head

Poisoning and caustic burns, or undesirable reactions caused by escaping hazardous materials

- Maintain the tightening torque of connections (see Hydraulic connections)
- Check tightening torque of head screws regularly
- > Only use pump if it is in perfect working order
- > Operate pump in accordance with the technical data

A DANGER

Danger of injuries and property damage caused by medium escaping from the discharge outlet

Poisoning and caustic burns, or undesirable reactions caused by escaping hazardous materials

- Place the pump in a suitable collecting vessel
- Never close the discharge outlet
- If medium escapes from the pump's discharge outlet, stop using the pump and contact your KNF dealer

Risk of fire from escaping flammable media

Escaping or leaking flammable media can produce a combustible environment.

- Take fire protection precautions that are appropriate for the transferred media
- Keep ignition sources away
- Keep suitable extinguishers accessible
- Use means of fire detection

WARNING

Danger of serious injuries or property damage caused by escaping media in the event of overpressure on the suction side

Excessive pressure on the suction side will result in uncontrolled escaping of the fluid, even if the pump is not operating

- Avoid overpressure on the suction side
- Fit a pressure control valve (see Section 6.3.3)









A CAUTION

Deviating metering accuracy

The metered volume may deviate from the set value due to gases, particles, or cavitation in the fluid system.

- Use a pre-filter, especially for media containing solid particles or fibers (see Chapter 13)
- Clean the valve disc/pump head (see Section 10.4)
- Ventilate the pump and system well
- Set the liquid type, for highly viscous and outgassing media, see Section 7.11

A CAUTION

Deviating metering accuracy

Transferring media containing particles larger than 70 µm may block the valves in the pump head.

- Use a pre-filter, especially for media containing solid particles or fibers (see Chapter 13)
- Monitor the metering
- Clean the valve disc/pump head (see Section 10.4)

2.6 Personal protective equipment

Special protective equipment must be used when performing certain tasks. These are indicated separately in the individual Chapters.

Where the wearing of protective equipment is concerned it is essential to observe the safety data sheet for the medium to be transferred.

The following symbols appear in the working area or in the operating instructions:

Safety gloves

to protect skin from contact with the transferred medium in accordance with the safety data sheet.



Safety glasses

to protect the eyes from contact with the transferred medium in accordance with the safety data sheet.

Face protection

to protect skin and eyes from contact with the transferred medium in accordance with the safety data sheet.







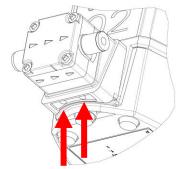


Fig. 1: Discharge outlet





2.7 Safety equipment

2.7.1 Safety diaphragm with intermediate flange

Danger of injuries and property damage caused by medium escaping from the discharge outlet

Poisoning and caustic burns, or undesirable reactions caused by escaping hazardous materials

- Place the pump in a suitable collecting vessel
- Check the discharge outlet regularly for leaks
- Never close the discharge outlet (see Fig. 1)
- If medium escapes from the pump's discharge outlet, stop using the pump and contact your KNF dealer

In the event that the pump diaphragm ruptures, the safety diaphragm prevents the transferred medium from penetrating into the pump. If the pump diaphragm ruptures, the transferred medium is channelled out of the pump through the discharge outlet (see Fig. 1).

2.7.2 Emergency Stop

Pulling out the mains plug acts as an Emergency Stop. If the pump is installed permanently, an Emergency Stop and a maintenance switch must be fitted.

2.7.3 Pressure control valve (optional)

In the event of overpressure on the suction side, the pump allows the medium to flow through it freely. The pressure control valve maintains a counterpressure of 0.5 bar and prevents the transferred medium from escaping. The pressure control valve is to be used on the pressure side (see Section 6.3.3)

2.8 Environmental protection considerations

All replacement parts should be properly stored and disposed of in accordance with the applicable environmental protection regulations. Ensure adherence to the pertinent national and international regulations. This especially applies to parts contaminated with toxic substances.

Dispose of any packaging material that is no longer required in an environmentally friendly manner. The packaging materials are recyclable.

Dispose of end-of-life equipment in an environmentally friendly manner. Use appropriate waste collection systems for the disposal of end-of-life equipment. Used pumps contain valuable recyclable materials.

2.9 Danger areas

When transferring corrosive, poisonous, ionising, inflammable or biological media, escaping medium may cause a hazard.

The pump danger area should be limited by safe workplaces. Safe workplaces should comply with the requirements of local work safety regulations.

2.10 Declaration of conformity

		K – Konformi K Declaration	
Hiermit erklärt der He Herewith the manufact			
KNF Flodos AG, Was	5	0 Sursee, Schw	/eiz
dass folgende Produ that the following prod			
Bezeichnung Denomination		Modell Model	Тур Туре
Labor Membran	-Dosierpumpen	FEM 1.02	.18S/. 18RC/. 18RCP
Labor Membran		FEM1.10	.18S2/. 18RC2/. 18RCP2
Labor Membran		UFEM 1.10	.18S2/. 18RC2/. 18RCP2
(Seriennummer siehe (Serial number see type			
comply with all relevan Richtlinie 2006/4 Directive 2006/42		llowing Directive: htlinie) tive)	/ Verordnungen entsprechen: s / Regulations:
ok negulation 5.1.		Machinery (Salet	у)
Richtlinie 2014/3 Directive 2014/30 UK Regulation S.I. Richtlinie 2011/6	0/EU (EMV) /EU (EMC) 2016/1091 Electroma 55/EU (RoHS)		
Richtlinie 2014/3 Directive 2014/30 UK Regulation S.I. Richtlinie 2011/6 Directive 2011/65	0/EU (EMV) /EU (EMC) 2016/1091 Electroma 5/EU (RoHS) /EU (RoHS) 2012/3032 Restriction rent	agnetic Compatibi n of the Use of Ce zugrunde geleg	lity ertain Hazardous Substances in Electrical and
Richtlinie 2014/3 Directive 2014/30 UK Regulation S.I. Richtlinie 2011/6 Directive 2011/65 UK Regulation S.I. Electronic Equipm Folgende harmonisie	0/EU (EMV) /EU (EMC) 2016/1091 Electroma 5/EU (RoHS) /EU (RoHS) 2012/3032 Restriction rent	agnetic Compatibi n of the Use of Ce zugrunde geleg	lity ertain Hazardous Substances in Electrical and
Richtlinie 2014/3 Directive 2014/30 UK Regulation S.I. Richtlinie 2011/6 Directive 2011/65 UK Regulation S.I. Electronic Equipm Folgende harmonisie The following harmoni	0/EU (EMV) /EU (EMC) 2016/1091 Electroma 55/EU (RoHS) /EU (RoHS) 2012/3032 Restriction ent ent ente Normen wurden ized standards have be Ausgabedatum	agnetic Compatibi n of the Use of Ce zugrunde geleg en applied: Titel Titel Title Dcchnische Dc	lity ertain Hazardous Substances in Electrical and It: ikumentation zur Beurteilung von Elektro- un
Richtlinie 2014/3 Directive 2014/30 UK Regulation S.I. Richtlinie 2011/6 Directive 2011/65 UK Regulation S.I. Electronic Equipm Folgende harmonisie The following harmoni Fundstelle Reference	0/EU (EMV) /EU (EMC) 2016/1091 Electroma 5/EU (RoHS) /EU (RoHS) 2012/3032 Restriction ent erte Normen wurden ized standards have be Ausgabedatum Release date	agnetic Compatibi n of the Use of Ce zugrunde geleg en applied: Titel Title Technische Dc Elektronikgeräter Sicherheit von	lity ertain Hazardous Substances in Electrical and It: humentation zur Beurteilung von Elektro- un n hinsichtlich der Beschränkung gefährlicher Stoffe Maschinen - Allgemeine Gestaltungsleitsätze
Richtlinie 2014/3 Directive 2014/30 UK Regulation S.I. Richtlinie 2011/6 Directive 2011/65 UK Regulation S.I. Electronic Equipm Folgende harmonisie The following harmoni Fundstelle Reference EN IEC 63000	0/EU (EMV) /EU (EMC) 2016/1091 Electroma 5/EU (RoHS) /EU (RoHS) 2012/3032 Restriction ent ent ent ent ent ent Ausgabedatum Release date 2018	agnetic Compatibi n of the Use of Ce zugrunde geleg en applied: Titel Title Technische Do Elektronikgeräter Sicherheit von Risikobeurteilung Elektrische Me	lity ertain Hazardous Substances in Electrical and It: humentation zur Beurteilung von Elektro- un n hinsichtlich der Beschränkung gefährlicher Stoffe Maschinen - Allgemeine Gestaltungsleitsätze g und Risikominderung ss-, Steuer-, Regel- und Laborgeräte - EMV
Richtlinie 2014/3 Directive 2014/30 UK Regulation S.I. Richtlinie 2011/65 UK Regulation S.I. Electronic Equipm Folgende harmonisie The following harmoni Fundstelle Reference EN IEC 63000 EN ISO 12100	0/EU (EMV) /EU (EMC) 2016/1091 Electroma 55/EU (RoHS) /EU (RoHS) 2012/3032 Restriction ent erte Normen wurden ized standards have be Ausgabedatum Release date 2018 2011	agnetic Compatibi n of the Use of Ce zugrunde geleg en applied: Titel Titel Technische Do Elektronikgeräter Sicherheit von Risikobeurteilung Elektrische Me Anforderungen - Sicherheitsbesti	lity ertain Hazardous Substances in Electrical and it: humentation zur Beurteilung von Elektro- un hinsichtlich der Beschränkung gefährlicher Stoffe Maschinen - Allgemeine Gestaltungsleitsätze und Risikominderung ss-, Steuer-, Regel- und Laborgeräte - EMV Teil 1: Allgemeine Anforderungen mmungen für elektrische Mess-, Steuer-, Regel- un
Richtlinie 2014/3 Directive 2014/30 UK Regulation S.I. Richtlinie 2011/6 Directive 2011/65 UK Regulation S.I. Electronic Equipm Folgende harmonisie The following harmoni Fundstelle Reference EN IEC 63000 EN ISO 12100 EN 61326-1 EN 61010-1 Bevollmächtigte Pers Authorised person to c	0/EU (EMV) /EU (EMC) 2016/1091 Electroma 5/EU (RoHS) /EU (RoHS) 2012/3032 Restriction ent ent ent Normen wurden ized standards have be Ausgabedatum Release date 2018 2011 2013 2010/A1:2019 son für die Zusamme ompile the relevant teo	agnetic Compatibi n of the Use of Ce zugrunde geleg en applied: Titel Titel Technische Do Elektronikgeräter Sicherheit von Risikobeurteilung Elektrische Me Anforderungen - Sicherheitsbestin Laborgeräte - Te	lity ertain Hazardous Substances in Electrical and t: humentation zur Beurteilung von Elektro- un hinsichtlich der Beschränkung gefährlicher Stoffe Maschinen - Allgemeine Gestaltungsleitsätze und Risikominderung ss-, Steuer-, Regel- und Laborgeräte - EMN Teil 1: Allgemeine Anforderungen mmungen für elektrische Mess-, Steuer-, Regel- un il 1: Allgemeine Anforderungen

3 Technical data

3.1 Items included in delivery

- SIMDOS metering pump
- Mains plug
- Operating manual
- Abbreviated instructions
- Connection kit: Hose and screw-in fittings
- Filter
- Remote control cable (RC and RCP version only)
- D-sub control cable (RCP version only)

3.2 Storage conditions

- Pumps must be stored in a dry place and protected from contamination.
- The storage temperature must be between 5°C and 40°C.
- The pump is supplied with protective caps. These protective caps must be fitted during storage.
- Store pump upright and with protection.

3.3 Pump head materials

The type designation **KT** stands for:

Assembly	Material ¹⁾	Medium contact
Pump head	PP	Yes
Valve plate / seals	FFKM	Yes
Diaphragm	PTFE-coated EPDM membrane	Yes
Included Filter	PEEK*	Yes
Housing	PP, TPE, PC	No

Tab. 1: KT materials

¹⁾ according to DIN ISO 1629 and 1043.1 *material for housing and filter mesh

The type designation **TT** stands for:

Assembly	Material ¹⁾	Medium contact
Pump head	PVDF	Yes
Valve plate / seals	FFKM	Yes
Diaphragm	PTFE-coated EPDM membrane	Yes
Included Filter	PVDF*	Yes
Housing	PP, TPE, PC	No

Tab. 2: TT materials

¹⁾ according to DIN ISO 1629 and 1043.1 *material for housing and filter mesh

The type	designation	FT	stands	for:
The type	acoignation		otunuo	101.

Assembly	Material ¹⁾	Medium contact
Pump head	PTFE	Yes
Valve plate / seals	FFKM	Yes
Diaphragm	FFKM	Yes
Included Filter	PEEK*	Yes
Housing	PP, TPE, PC	No

Tab. 3: FT materials

¹⁾ according to DIN ISO 1629 and 1043.1 *material for housing and filter mesh


The type designation ST stands for:

Assembly	Material	Medium contact
Pump head	1.4404	Yes
Valve plate / seals	FFKM ¹⁾	Yes
Diaphragm	PTFE-coated EPDM membrane	Yes
Included Filter	PEEK*	Yes
Housing	PP, TPE, PC ¹⁾	No

Tab. 4: ST materials


¹⁾ according to DIN ISO 1629 and 1043.1 *material for housing and filter mesh

3.4 Product key

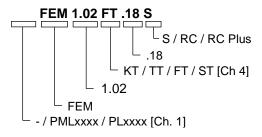

Description	Meaning	
PML / PL	Customer-specific designs	
FEM	Flodos Electronic Metering (designation for metering pumps)	
1.02	Pump type	
KT / TT / FT / ST	Head material	
.18	Laboratory device with housing	
S / RC / RC Plus	Swithout remote controlRCwith remote control (Remote Control)RC Pluswith RS 232 interface	

Tab. 5: Product key

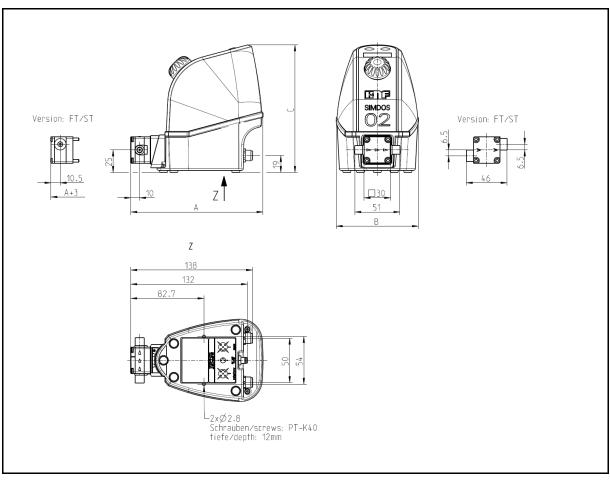

3.5 Dimensions

Fig. 2: Mounting measurements (measurement tolerances as per DIN ISO 2768-1, tolerance class V)

Parameter	Value (metric)	Value (imperial)
Width (A)	150 mm	6.102 inches
Depth (B)	93 mm	3.740 inches
Height (C)	144 mm	5.708 inches
Weight of pump ¹	0.9 kg	1.98 lb

Tab. 6: Dimensions

Type plate

The type plate is on the bottom of the device. It contains the following information:

- Manufacturer's address
- Type
- Serial Number
- Design pressure
- CE and UKCA mark
- Year of manufacture
- Electrical connection data

¹ Differences are possible depending on the project

3.6 Installation location requirements

The pump is designed to be operated in an upright position.

NOTE An unsuitable location may cause damage to the device \triangleright Install in a dry place Protect the installation location of the pump from water in the form of rain, spray, splashes and drips > Do not operate the pump in an aggressive or inflammable environment Choose a safe location (a level, flat surface) for the pump \geq Place the pump in a suitable collecting pan > If the stability of the pump cannot be ensured, fasten the pump in place using two screws > Make sure that the hose connections can be fitted without strain on the hoses > Protect the pump against shocks, impacts and strong vibrations

3.7 Electrical connections and performance data

3.7.1 Power supply unit

Value
100 – 240 V AC +/- 10 %
50 - 60
40
24
IP 40
Electronic overload protection

Tab. 7: Electrical data power supply unit

3.7.2 Pump

Parameter	Value
Pump DC voltage [V]	24
Max. current consumption, DC RMS 24 V [A]	0.4
Max. short-term peak current [A]	1.7
Pump protection	Electronic overload protection
Pump protection type	IP 65

Tab. 8: Electrical data pump

Parameter	Value
Analog input	
Signal range, (100%) for 1 – 100% of full scale	0 – 5 V, 0 – 10 V, 0 – 20 mA, 4 – 20 mA
Signal range, (30%) for 0.3 – 30% of full scale	0 – 5 V, 0 – 10 V, 0 – 20 mA, 4 – 20 mA
Signal range, (15%) for 0.15 – 15% of full scale	0 – 5 V, 0 – 10 V, 0 – 20 mA, 4 – 20 mA
Input resistance [Ω]	13.9 kΩ at 0 – 10 V 13.9 kΩ at 0 – 5 V 470 Ω at 4 – 20 mA 470 Ω at 0 – 20 mA
Electric strength [V] TTL	24 V DC
Digital input	
Signal range	Pull up at 24 V
Voltage [V] TTL	24 V DC
Low level (ON)	< 0.8 V = low
High level (OFF)	> 2.0 V = high
Digital output	
Voltage open collector [V] TTL	24 V DC
Load capability open collector [mA] TTL	10 mA

3.8 Remote control (RC version only)

Tab. 9: Remote control

3.9 Remote control via RS 232 (RC Plus version only)

Parameter	Value
RS 232 interface	
TX – RS232	
RX – RS232	
Ground	
Settings (factory)	
Baud rate	9600 Kbit/s
Data	Bits 8
Parity	No
Stop	Bit 1
Mode	Half duplex

Tab. 10: RS 232 interface

3.10	Flow	rates
------	------	-------

Parameter	Value
Flow rate [ml/min]	0.03 - 20
Max. metering volume [ml]	0.03 - 1000
Metering time [mm:ss] [hh:mm]	1s – 99h 59min
Permissible pressure [bar g]	6
Suction head [mWG]	2

Tab. 11: Hydraulic ratings

NOTE

The pump has been calibrated to its nominal flow rate in the factory. The specified flow rate is for water at 20°C with free discharge flow. The flow rate may differ from the calibrated value depending on the medium, its viscosity, density and temperature, pressure and installation type (e.g. cross-section constrictions).

 For accurate metering, calibrate the pump in the system (see Section 7.12)

3.11 Other parameters

Parameter	Value
Permissible ambient temperature [°C]	+5 to +40
Service life [h]	10,000 (100 million strokes) or 5 years
Noise level	< 40 dBA

Tab. 12: Other parameters

3.12 Hydraulic connections

Pump type	Connection type
FEM 1.02 KT.18	Internal thread UNF 1/4" - 28
FEM 1.02 TT.18	Internal thread UNF 1/4" - 28
FEM 1.02 FT.18	Internal thread UNF 1/4" - 28
FEM 1.02 ST.18	Internal thread UNF 1/4" – 28

Tab. 13: Hydraulic connections

3.13 Transferred medium

Parameter	Value
Permissible medium viscosity [cSt] ²	150
Ideally free of solids, max. particle size [µm]	< 70
Permissible media temperature ³ [°C]	+5 to +80

Poisoning and caustic burns

If corrosive, poisonous, ionising or biological media come into contact with skin, eyes and the respiratory tract they may cause serious injuries

- Observe the safety data sheets for the media to be transferred
- > Use personal protective equipment
- > Operate the pump in a suitably protected area
- Check the resistance of any parts which come into contact with the fluid (see chemical resistance list or consult your local KNF dealer)

Damage caused by ionising radiation

Radioactive materials cause serious damage to health and reduce the service life of the pump.

- Decontaminate the pump
- Wear personal protective equipment
- Mark the pump

NOTE

If it cannot be ensured that the transferred medium is free of solids, $a < 70 \mu m$ filter with a sufficiently large filter area must be fitted upstream of the pump to prevent it from being damaged (see Chapter 13).

NOTE

Before using a medium, check the compatibility of the materials of the pump head, pump housing, diaphragm and valves with the medium.

- Check resistance, see chemical resistance list and Section 4.3
- Clarify resistance with KNF





² Using the fluid type setting "Visc500cSt" allows media with a viscosity of 500 cSt to be transferred at a limited flow rate. (see Section 7.11)

³ Applies to water and low-reactivity media. For aggressive media, resistance is to be checked at a higher medium temperature.

3.14 Accuracy/ reproducibility

The accuracy of the metering pump is appropriately characterised by the maximum absolute error A and the maximum relative error B:

➔ max. inaccurac	$y \le A + B$
------------------	---------------

Error	Value
Absolute error A	≤ ± 1‰ nominal value
Relative error B	≤ ±1.9% of setting value
Max. deviation	≤ A+B
Mean accuracy	+/- 2%
Mean reproducibility ⁴	+/- 1%

Tab. 14: Error A and error B

Example: SIMDOS 02, nominal value of 20ml/min

For a setting of 5ml/min, the accuracy should therefore be within the following error limits:

Deviation

- \leq (±1‰ x 20ml/min) + (±1.9% x 5ml/min) \leq (±0.02ml/min) + (±0.095ml/min)
 - = (±0.02mi/mir) · (±0.0
 - ≤ ±0.115ml/min
 - ≈ ±2% of setting value

A CAUTION

Deviating metering accuracy

The metered volume may deviate from the set value due to gases, particles, or cavitation in the fluid system.

- Use a pre-filter, especially for media containing solid particles or fibers (see Chapter 13)
- Clean the valve disc/pump head (see Section 10.4)
- Ventilate the pump and system well
- Set the liquid type, for highly viscous and outgassing media, see Section 7.11

A CAUTION

Deviating metering accuracy

Transferring media containing particles larger than 70 µm may block the valves in the pump head.

- Use a pre-filter, especially for media containing solid particles or fibers (see Chapter 13)
- Monitor the metering
- Clean the valve disc/pump head (see Section 10.4)



⁴ Under constant environmental conditions and with the same hose configuration

NOTE

Greater inaccuracies may occur for fluids with a viscosity of > 150 cSt or with a tendency to gassing out

- Set "fluid type" (see Section 7.11)
- Calibrate the pump in the system (see Section 7.12)

NOTE

The flow rate may differ from the calibrated value for some applications, according to the medium, its viscosity, density and temperature, pressure and installation type (e.g. cross-section constrictions)

 For accurate metering, calibrate the metering pump in the system (see Section 7.12)

4 Assembly and function

4.1 Metering pump structure

- 1 STOP button
- 2 START button
- 3 Display
- 4 Control knob
- 5 Inlet
- 6 Pump head
- 7 Outlet
- 8 Power supply connector
- 9 Remote Control connector (RC version only)
- 10 Connector for serial communication (RC Plus version only)

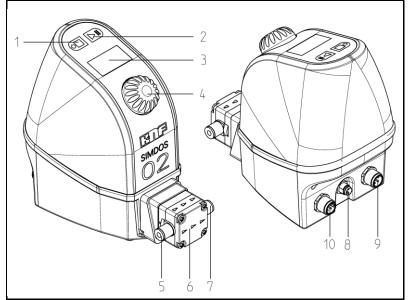


Fig. 3: Diaphragm pump

4.2 Operating principle

- 1 Outlet valve
- 2 Inlet valve
- 3 Working chamber
- 4 Diaphragm
- 5 Eccentric
- 6 Connecting rod
- 7 Pump drive

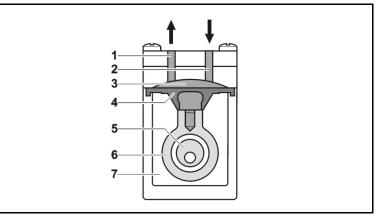


Fig. 4: Pump assembly

Principle:

The diaphragm liquid pumps are based on reciprocating displacement pump technology. The elastic diaphragm (4) is moved up and down by the eccentric (5) and the connecting rod (6). In the downward stroke it aspirates the medium to be transferred via the inlet valve (2). In the upward stroke, the diaphragm pushes the medium out of the pump head via the outlet valve (1). The diaphragm hermetically seals off the working chamber (3) from the pump drive (7). Patented drive system

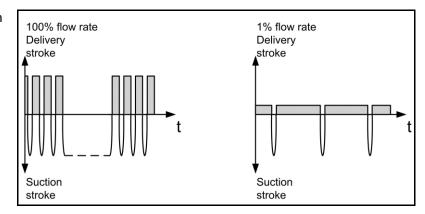



Fig. 5: Low-pulsation operating principle

SIMDOS[®] diaphragm metering pumps are fitted with a patented drive system. The suction stroke is always driven at maximum speed, but the exhaust process is varied according to the selected speed, ensuring maximum possible uniformity of discharge during the entire stroke (see Fig. 4). This results in a low-pulsation, quasicontinuous transfer process. This is the key difference between SIMDOS[®] metering pumps and diaphragm metering pumps with linear magnet diaphragm drive or with conventional eccentric diaphragm drive.

The characteristic curve for the suction and exhaust strokes can also be adapted to the medium by setting the appropriate fluid type (see Section 7.11).

4.3 Valves

The SIMDOS is supplied with DuPont[™]Kalrez^{®5} valve disks as standard, but can also be supplied with valve disks made of CHEMRAZ^{®6} as an option.

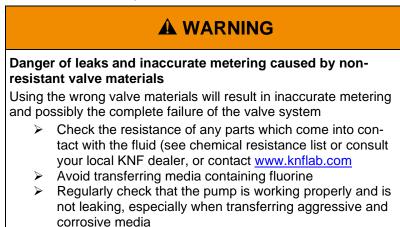



⁵ DuPont™Kalrez® is a registered trademark of E.I. du Pont de Nemours and Company or one of its subsidiaries. All rights reserved.
⁶ CHEMRAZ[®] is a registered trademark of Greene, Tweed & Co. All rights reserved.

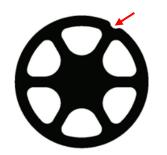


Fig. 6: Kalrez[®], standard valve

4.3.1 Standard valve, DuPont[™]Kalrez[®]

The SIMDOS is supplied with DuPont™Kalrez[®] valve disks as standard. This material is resistant to most chemicals.

The optional CHEMRAZ[®] Valve Kit (see Section 13.1) is required for the liquids in Tab. 15.

- The resistance of the individual materials can be found in the chemical resistance list provided.
- > In case of doubt, please contact your local KNF dealer.

NOTE

The valve material can be identified from the geometry of the valve disk.

- DuPont™Kalrez[®] has one groove in its outer contour (see Fig. 6)
- CHEMRAZ[®] has six grooves in its outer contour (see Fig. 7)
- If CHEMRAZ[®] valves are used, the pump is to be labelled accordingly
- The instructions in Section 10.4 are to be followed when removing the valves

4.3.2 CHEMRAZ[®] Valve Kit

Experience shows that $\mathsf{CHEMRAZ}^{\texttt{®}}$ values should be used for the following liquids:

Solvents		
Tetrahydrofuran (THF)	C ₄ H ₈ O	
Dimethyl sulfoxide (DMSO)	C ₂ H ₆ OS	
Dimethyl formamide (DMF)	C ₃ H ₇ NO	
Methyl ethyl ketone (MEK)	C ₄ H ₈ O	
Cyclohexane	C ₆ H ₁₂	
Concentrated acids (75%)		
Hydrofluoric acid	HF	
Nitric acid	HNO ₃	
Sulphuryl dichloride	SO ₂ CL ₂	
Trifluouroacetic acid	C ₂ HF ₃ O ₂	

Tab. 15: Excerpt from the chemical resistance list for the CHEMRAZ[®] Valve Kit

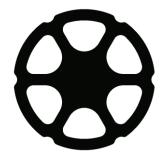


Fig. 7 Chemraz[®] valve, optional

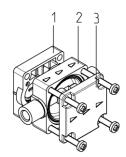


Fig. 8: Pump head


4.4 Pump diaphragm

Operating controls

The pump diaphragm is covered with a layer of PTFE to ensure good chemical resistance. Type "FT" uses a pure FFKM diaphragm.

4.5 Pump head

The pump head comprises an intermediate plate (1), a connecting plate (2) and a head plate (3). The intermediate and connecting plates come into contact with the liquid, so the resistance of the corresponding head materials must be established.

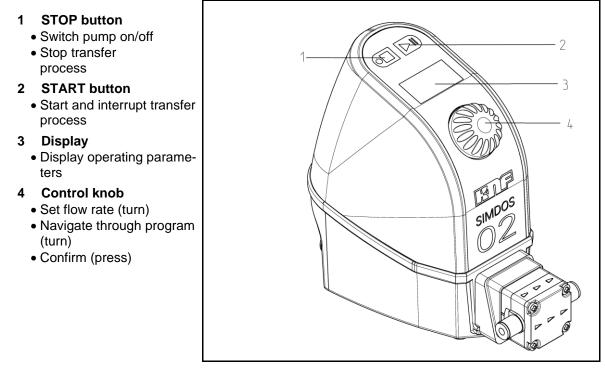

4.6

Fig. 9: Operating controls

	The main display shows the principal pump settings.	
 Specified flow Cycle metering mode Current cycle Programmed cycles Autostart Remote control Analog input Digital input 1 Digital input 2 Operating mode display Flow setting unit Volumeter Time counter 	$\begin{array}{c ccccccccccccccccccccccccccccccccccc$	
Specified flow and unit	<i>Fig. 10: Main display</i> Displays the flow rate set (1) if the unit (11) has been set to "ml/min". Displays the metering volume (1) if the unit (11) has been set to "ml".	
Time counter	 Displays the pump running time. Downward count (13) if a timeout (metering) has been programmed. Time remaining display. 	
Volumeter	Display of delivery volume (12) since last pump start-up.	
Remote control	Displays whether pump external drive has been activated, and the external drive condition (6, 7, 8, 9).	
Autostart	Displays whether pump has been set to Autostart (5).	
Operating condition	Displays whether the pump is operating, stopped or paused (10).	
Cycle metering mode	If this display (2) is active, the programmed	
	number of volumes (3) is metered. During operation the current cycle number (3) is incremented until the programmed cycle number (4) is reached.	
	 During metering, the transfer time remaining for the current dose is displayed. 	
	 During the pause the pump counts down the pause time remaining. 	

4.6.1 Display

The main display shows the principal pump settings.

Display	Function
20.0 ml/min minsec 0.0 ml	Flow rate 0.03 – 20 ml/min Metering volume 0.03 – 1000 ml and metering time timeout
Prime	High-speed operation for priming or emptying
Fluid type S	Set the pump for the characteristics of the medium to be trans- ferred
	S = Standard: aqueous media
	D = Volatile: media with a low boiling point
	V = Visc100 cSt: media up to 100 cSt
	H = Visc500 cSt: media up to 500 cSt
Cal. 100%	To calibrate the pump
Cycle Off	Set cycle metering mode (several volumes) Quantity: Repetitions (Off, 2999, Inf) Pause: pause time (1s99 min 59s)
System →	Switch to system menu


4.6.2 Main menu


Display	Function
Back ←	Switch to main menu
English	select language
A_in Aı	Analog input of pump* ■ Off ■ 0 – 5 V, 0-10 V, 0 – 20 mA, 4 – 20 mA
Range 100	 % Flow rate range selection for analog input* 100% 1 – 100% 0.20 – 20 ml/min 30% 0.3 – 30% 0.06 – 6 ml/min 15% 0.15-15% 0.03 – 3 ml/min
ln1 c	 Digital input 1:* Off Start/Stop using level or pulse signal
ln2 c	 Digital input 2:* Off Reset or high-speed operation
Out Alar	 m Digital output:* Error/alarm End of metering volume Volume pulses Motor pulses
Autostart c	 Set autostart Off On (pump will start running as soon as mains voltage is applied)
Contrast	40 Set display contrast
Reset Ne	in Reset pump to factory settings
Maintenanc.	Maintenance position for diaphragm replacing
Back	Switch to main menu
1	ΝΟΤΕ
	* RC functions are available for RC pumps only (see Section 3.4). See Chapter 8 for further details of the external drive system.

4.6.3 System menu

4.7 Operating modes

Continuous transferring In this case, the device transfers at a constant flow rate.

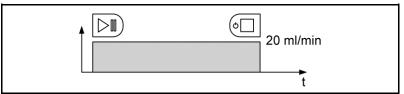

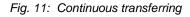

Continuous transferring with timeout

In this case, the device transfers at a constant flow rate, then switches off after a specified time.

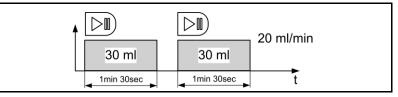


Fig. 12: Continuous transferring with timeout

Metered volume

Transferring of a specified volume in a specified time.

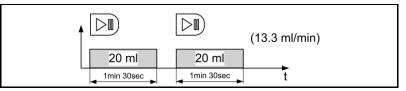


Fig. 13: Metered volume with timeout

Cycle metering mode metering several volumes

The pump meters n repetitions of a defined metering cycle with pause time.

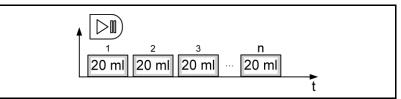


Fig. 14: Cycle metering mode

5 Shipment

5.1 Checking delivery

Prior requirements

 Device delivered and packaged

Procedure

- ure 1. Check delivery for visible damage
 - 2. Check that delivery is correct and complete using the packing list
 - 3. In the event of a discrepancy, register a complaint.
 - 4. Keep original packaging in case you have to return the item

5.2 Registering a complaint

Prior requirements Device damaged or incomplete

Procedure

- 1. If there are any visible signs of damage, only accept the delivery with reservations.
 - 2. Damage and complaints of any kind should be recorded on the delivery note
 - 3. Notify the manufacturer of any damage and complaints

5.3 Arranging for return

Prior requirements • Device damaged or incomplete

Procedure

- 1. Contact KNF before returning the device
- 2. Flush and decontaminate pump
- 3. Fill out decontamination declaration and attach to the pump
- 4. Send the pump in its original packaging to a service location or one of our product centres
- 5. The pump can be sent by parcel service or post in its original packaging

A DANGER

Danger of poisoning, chemical burns, or radioactive contamination through contact with pump parts that are not decontaminated

Contact with pump parts that are not properly decontaminated will result in poisoning, chemical burns, or radioactive contamination.

- > Use personal protective equipment
- Flush and decontaminate the pump until the possibility of danger can be eliminated.
- Label any pumps that are used with particularly dangerous media
- Only decontaminated pumps may be sent



NOTICE

KNF retains the right to refuse repair of pumps that are used to move hazardous materials.

This applies particularly to the following materials:

- Poisons
- Radioactive media
- Elementary halogens (e.g. bromine)

6 Installation and initial start-up

6.1 Safety

A WARNING

Danger as a result of incorrect installation

Injuries or damage to equipment caused by escaping media

- Follow installation instructions
- > After assembly, test the system with a harmless liquid
- > Only use pump if it is in perfect working order

Danger of serious injuries or property damage caused by escaping media in the event of rupturing of the pump diaphragm

Poisoning and caustic burns caused by escaping media as a result of rupturing of the pump diaphragm. If the pump diaphragm ruptures, the medium is channelled out of the pump through the discharge outlet.

- When transferring critical media, operate the pump in a collecting tank
- Check discharge outlet for leaks

Danger of serious injuries or property damage caused by escaping media in the event of overpressure on the suction side

Overpressure on the suction side will result in uncontrolled escaping of the fluid, even if the pump is not operating.

- Avoid overpressure on the suction side
- Fit a pressure control valve (see Section 6.3.3)

The following requirements apply to all the activities described in this chapter:

Prior requirements

- Device is in perfect working order
- The system is operated with water
- Must be assembled by specialist personnel

Personal protective equipment

Not required

Personnel







6.2 Installation

Before installation, store the pump at the installation location to bring it up to ambient temperature.

NOTE

For problem-free, safe operation the following instructions should be followed:

- Install in a dry place.
- Protect the installation location of the pump from water in the form of rain, spray, splashes and drips.
- > Do not operate the pump in an aggressive or inflammable environment.
- > Choose a safe location (a level, flat surface) for the pump.
- If the stability of the pump cannot be ensured, fasten the pump in place using two screws.
- Make sure that the hose connections can be fitted without strain on the hoses.
- Protect the pump against shocks, impacts and strong vibrations.

6.2.1 Connect mains plug

Risk of electric shock

- Only operate the pump with the power supply unit supplied
- Connect the cable from the mains plug to the pump's power supply socket
- Connect the mains plug into a correctly installed and properly earthed mains socket
- Ensure that the mains plug is protected from water splashes
- > Do not open pump housing or power supply unit
- Disconnect pump from power supply during maintenance work
- > If pump is installed permanently, fit a maintenance switch

When connecting the device to a power source, the relevant norms, directives, regulations and technical standards must be observed.

- 1. Use suitable adapters in the power supply unit
- 2. Connect power supply unit to pump

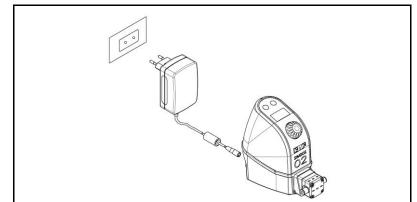


Fig. 15: Electrical connection



6.2.2 Connect filter



^ **_** _ _ _ _ _ _ _ _ _ _

Danger of excessive banking-up pressure (resistance) in the filter as a result of high particle concentration in the medium or a mesh size which is too small

This reduces the flow rate and causes errors in metering accuracy

- Use filters with a sufficiently large mesh size
- Filter fluid beforehand where appropriate
- Check filter penetration regularly, cleaning or replacing the filter where necessary (see filter data sheet)

Danger of blocked filters as a result of excessive particle concentration in the medium

This reduces the flow rate and causes errors in metering accuracy

- Filter fluid beforehand where necessary
- Check filter penetration regularly, cleaning or replacing the filter where necessary (see filter data sheet)

Many media contain solids or fibres that are not visible to the naked eye. In order to maintain reliable pump operation, a filter is to be used on the suction side.

- 1. Connect the filters supplied on the suction side (for Order No. see Section 13.4)
- 2. Set up the system in the following steps
- 3. Rotate pump 90° anticlockwise
- 4. Prime pump so that the air can escape from the filter.
- 5. If there are no air bubbles in the system, the pump can be put into operation.

NOTE

Air in the pump head causes inaccurate metering.

- 1. Empty pump
- 2. Refill pump

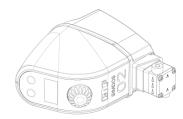


Fig. 16: Prime pump vertically

6.2.3 **Connect hydraulics**

Connected

1.

- components
- handle the hydraulic characteristics of the pump (see Chapter 3). Hoses
 - 2. Only use hoses that are suitable for the maximum operating pressure of the pump (see Section 3.10).
 - Only use hoses that are sufficiently chemically resistant to the 3. liquids being transferred.

Only connect components to the pump that are designed to


- Customer-specific pumps (PL, PML)
- The connections described below apply to standard products. 4. Different connections may apply for customer-specific projects (PML or PL).

NOTE

Arrows on the pump head indicate the flow direction.

FEM 1.02 - Screw-in connection fitting with internal thread

- Connector 1
- 2 Clamp ring
- Hose connector 3
- Hose 4

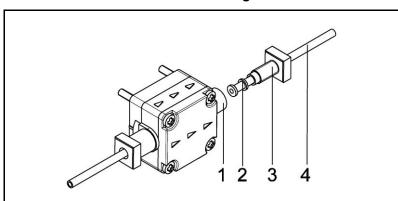

- Fig. 17: Hose connection UNF 1/4"-28
- 1. Remove protective caps
- 2. Push clamp ring (2) onto flanged hose
- 3. Push hose connector (3) onto flanged hose
- 4. Screw hose connector (3) into connector (1)
- 5. Hand-tighten hose connector (3), tightening a little more if necessary
- 6. Connect the suction and pressure lines. (Recommendation: UNF 1/4"-28 flanged connection with 1/8" hose, see accessories list in Section 13.2)
- 7. Check that the hoses (1) and transition joints (hose connector/hose) are connected correctly and securely
- 8. Check that the system is leak-tight (see Section 10.5)

6.3 System setup

6.3.1 Recommended setup

In order to maintain reliable pump operation, the following setups should be considered.

Description of symbols for the following diagrams:

Symbol	Meaning	Symbol	Meaning
\bigcirc	Pump	\mathbb{X}	Pressure control valve (adjustable)
\diamondsuit	Filter		Storage container

Tab. 16: Description of symbols

6.3.2 The storage container is level with the pump or underneath it

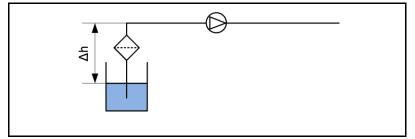


Fig. 18: Storage container underneath pump

In order to achieve the specified metering accuracy (see Section 10.5) the following instructions should be followed:

- Select suction head lower than Δh=1 mWG
- Keep suction hose as short as possible
- Select a suction hose with a large cross-section
- Vent the pump
- For constant results, keep the level of liquid constant (use a container with a large base)



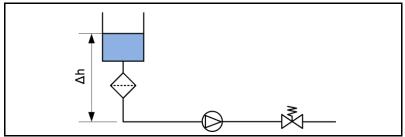


Fig. 19: Storage container above the pump.

Overpressure on the suction side reduces the accuracy of the pump

In order to avoid uncontrolled escaping of the medium, a pressure control valve must be used on the pressure side (see Fig. 19).

NOTE

- > Set up the system as shown in Fig. 19
- Suitable pressure control valves are available as accessories (see Section 13.3)
- Set the pressure control valve to 0.5 bar. This allows a Δh of less than 4.5 m to be compensated for

6.3.4 Metering volatile media

For precise metering of liquids with a low vapour pressure, such as solvents, the system should be set up as shown in Fig. 20

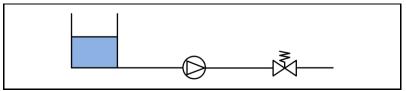


Fig. 20: Volatile media

- 1. Set up the system as shown in Fig. 20
- 2. The pressure control valve should be set to 0.5 bar
- To avoid cavitation, set the fluid type "Volatile" or "Visc500cSt" (see Section 7.11)

6.3.5 Protection against overpressure

In order to protect the system against overpressure, a safety valve must be fitted immediately downstream of the pump (see Fig. 21).

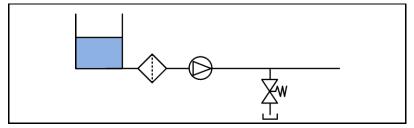


Fig. 21: Overpressure protection



6.4 Initial start-up

Danger of bursting of the fluid system as a consequence of overpressure

The pump builds up pressure. In a closed system the max. permitted operating pressure may be exceeded, thus causing damage to the pump and/or the system

- In the case of parts that are in contact with the fluid, only select those that are designed for the pump's operating pressure as a minimum
- Do not transfer against shut-off devices or closed systems, or
- install pressure relief/safety valves and set them to 6 bar (in the KNF product range)

Danger of unexpected chemical reactions with water

Water residues left in the pump from testing in the factory may react with the transferred medium.

 Before putting the pump into service, flush it with a noncritical medium

Before switching on the pump, check the following points:

Prerequisites for start-up

- All hoses attached properly
- Specifications of the power supply must correspond with the data on the pump and mains plug type plates.
- The pump outlet must be clear of any obstruction.
- All cables properly connected

Tab. 17: Prerequisites for start-up

- 1. Only operate the pump under the operating parameters and conditions described in Section 3.10.
- 2. Make sure the pump is used properly (see Section 2.1).
- 3. Avoid improper use of the pump (see Section 2.2).
- 4. Observe all safety precautions (see Chapter 2).

6.4.1 Test the system

In order to prevent damage, the safety, leak-tightness and functioning of the test set-up should be checked using a suitable harmless medium, such as water.



7.1 Safety

Danger of injuries and property damage caused by hazardous materials

Poisoning and caustic burns, or undesirable reactions caused by escaping hazardous materials

- Observe the safety data sheets for the media to be transferred
- > Determine the resistance of the head materials
- Maintain the tightening torque of connections (see 6.2)
- Check tightening torque of head screws regularly (see Section 10.4)
- Check that pump and system are leak-tight
- Service pump regularly
- Only use pump if it is in perfect working order
- > Operate pump in accordance with the technical data

A WARNING

Danger of undetected rupturing of diaphragm

The diaphragm may be damaged by aggressive media, and the medium may escape from the discharge outlet

- > Operate the pump in a collecting pan
- Flush pump thoroughly after use or if it is not to be used for an extended period (see Chapter 10.2)
- > Check the discharge outlet regularly for leaks
- Observe the service lives of diaphragms and other elastomer parts (see Maintenance)

A WARNING

Danger of splashing media

If corrosive, poisonous, ionising or biological media come into contact with skin, eyes and the respiratory tract they may cause serious injuries

- Observe the safety data sheets for the media to be transferred
- Check that the system is leak-tight before use
- > Use personal protective equipment

A CAUTION


Danger of cavitation or particles in the fluid circuit

Errors in metering accuracy

- > Use a prefilter for media containing particulate matter
- Clean valve disk/pump head (see Section 10.4)
- Vent the pump and system thoroughly
- Set the type of liquid in the case of high-viscosity and volatile media (see Section 7.11)





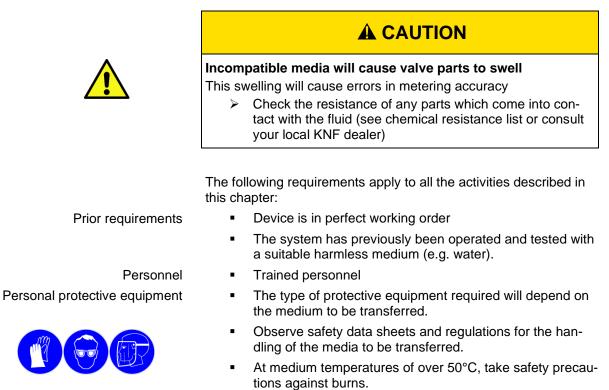

- When transferring aggressive, biological, inflammable or
- ionising media, a suitable work station with a safety and extraction system is to be used.

7.2 Switching pump on / off

The default setting is for the pump to remain off when connected to the mains power supply.

Switch on

1. Hold the "STOP" button down for at least 2 seconds

- ➔ the pump will now switch itself on
- → The display lights up, and the pump is ready for use.

Switch off

- 1. Hold the "STOP" button down for at least **2 seconds**.
 - → The pump will now switch off completely.

NOTE

When the pump is stopped, after ten minutes without any user input it switches into standby condition (display goes dark). However, the pump is still ready for use immediately when required.

- Press "STOP" or any other key to reactivate the pump from standby condition.
- The external drive reactivates the pump from standby as soon as a control signal is present.

L.

Special tools required

Fig. 22: Switching on the pump



Fig. 23: Switching off the pump

7.3 Entering settings

> Turn the control knob to scroll the display up or down.

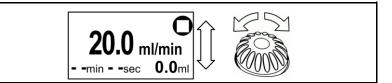


Fig. 24: Moving the operating menu up / down

The lines in the middle of the display are enclosed in a frame, and can now be selected with the control knob.


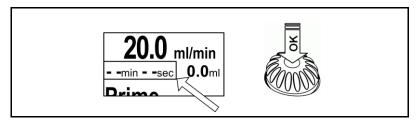


Fig. 25: Select the framed lines by pressing the control knob

The selected settings are marked with an inverted display. Now turn the control knob to change the value as required.

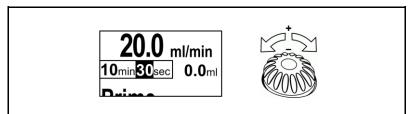

- Fig. 26: Inverted display of editable values
 - Press on the control knob to confirm your input, and move to the next value in the same line, if any.

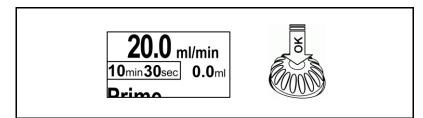


Fig. 27: Confirming the setting value by pressing the control knob **Ending the setting operation:**

Method 1

- 1. Keep pressing the control knob until no values are selected (values marked with inverted display).
- 2. Turn the control knob until the main display can be seen at the top of the operating menu.
- 3. Press the "STOP" button:
 - ➔ The pump will stop when this operation is carried out. This ends the input operation, and the display switches back to the main display.



Fig. 28: Specified flow

20 0	
20.0 m	nl/min
	0.0
minsec	0.0 ml

Fig. 29: Flow unit

Method 2

- 1. Press the "START" button: this ends the input operation, and the display switches back to the main display.
 - → The pump starts when this operation is carried out.

7.4 Setting the operating mode

- 1. Press control knob, Fig. 28 appears
 - ➔ The transfer value can be set by turning the control knob
- 2. Press the control knob to confirm the input
- 3. Turn control knob
- 4. Press control knob, Fig. 29 appears
 - ➔ The following operating modes can be set by turning the control knob.
 - [ml/min] for transferring at a fixed flow rate
 - [ml] for metering a volume
- 5. Press the control knob to confirm the input

Input	Unit	Setting range
Flow rate	Millilitres per minute[ml/min]	0.03 - 20.0
Metering volume	Millilitres [ml]	0.03 - 1000.0

Tab. 18: Operating mode setting ranges



Fig. 30: Prime

7.5 Priming the pump

This function is for rapidly priming and emptying the pump head and hoses.

- 1. Press the control knob and hold in place for the duration of the priming / emptying operation.
 - The pump will run at a higher speed during this time.

IMPORTANT

The priming function interrupts any metering operation that is currently in progress.

NOTE

Air bubbles in the hose system prevent precise metering.

- > Prime the system until no more bubbles are visible.
- > Avoid bubble formation in the fluid
- Hold pump with filter in vertical direction of flow and prime (see Section 6.2.2)

7.6 Starting the pump

- 1. Press the "START" button.
 - → The unit will begin transferring. The triangle symbol will appear in the display.

NOTE

An externally actuated pump will only start if the external control signals allow a pump start-up (see Chapter 8).

7.7 Interrupting a transferring operation

- 1. Press the "START" button
 - ➔ The transferring operation is interrupted. The pause symbol will appear in the display.

NOTE

In "Pause" operating mode:

- > when restarted, any interrupted transfer is continued.
- > the pump responds to external control signals.



Fig. 31: Starting the pump

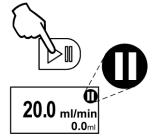


Fig. 32: Interrupting a transferring operation

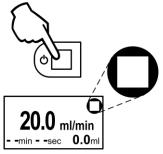


Fig. 33: Stopping the transferring operation

Master Stop

7.8 Stopping the transferring operation

- 1. Press the "STOP" button,
 - → The transferring operation is interrupted. The pause symbol will appear in the display.
- Press the "STOP" button again 2.
 - → The pump will now switch to "Stop" condition, and reset the time counter and volumeter.

NOTE

The first time the user presses the "STOP" button puts the working pump into Pause condition.

The pump goes into "Stop" condition only when the "STOP" button is pressed for the second time.

Stopping the pump with the "STOP" button is called a "Master Stop".

NOTE

Transferring after a Master Stop is only possible if:

- the Master Stop is manually cancelled using the "START" button:
- or if a "Reset" command is present at digital input 2.

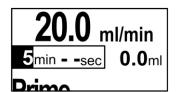


Fig. 34: Time setting in seconds

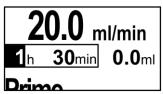


Fig. 35: Time setting in hours

7.9 Transferring fluid

- 1. Specify the required transfer value (see Section 7.4)
- 2. Set operating mode to [ml/min] (see Section 7.4)
- 3. Press the control knob to confirm the input
- 4. Press the "START" button
 - → Pump transfers at the desired flow rate

NOTE

When operating:

- a timeout can be programmed.
 - ➔ The pump transfers at a set flow rate until the time has elapsed
- if a timeout is not specified (see Fig. 29),
 - ➔ the pump transfers at the set flow rate until stopped by a Pause or Stop command
- the time counter can be switched to hours and minutes. To do this, set the minute setting to more than 59 min

Timeout with time counter

The timeout function can be configured as follows.

Input	Meaning	Setting range
min sec	Timeout not active	
mm min ss sec	Timeout active	1sec – 99h 59min

Tab. 19 Timeout

7.10 Setting metering operation

- 1. Specify the required transfer value (see Section 7.4)
- 2. Set operating mode to [ml] (see Section 7.4)
- 3. Press the control knob to confirm the input
- 4. Press the "START" button
 - ➔ The pump transfers the desired volume

NOTE

When metering a volume [ml]:

- the permissible setting range for the time counter is restricted so that the flow rate will not exceed or fall below the permissible flow rate for the pump.
- the time counter cannot be deactivated. The setting [- min - - sec] is not possible.

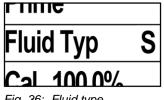


Fig. 36: Fluid type

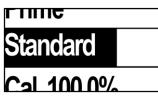


Fig. 37: "Standard" fluid type







7.11 Setting the fluid type

The "fluid type" setting is used to adjust the pump characteristic curve for different types of fluid.

- Turn control knob until "Fluid Type" appears in the selection 1. window
- 2. Press control knob
 - → The current setting is shown by character on the righthand edge

The following settings are available for selection:

Fluid type: Standard

Symbol: S

This is the optimum setting for all media with a viscosity similar to that of water. A uniform flow characteristic curve is obtained with short suction times and appropriate linearisation.

Fluid type: Volatile

Symbol: D

This is the appropriate setting for media with a low boiling point. As these media are sensitive to any pressure drop and deceleration phenomena:

- The pump suction rate is reduced
- The transferring characteristic curve reduces fluid acceleration and deceleration

Fluid type: Viscosity up to 100 cSt

Symbol: V

The setting "Visc100cSt" is for media with a viscosity of up to approx. 100 cSt. This setting alters the characteristics of the pump as follows:

- Suction rate is reduced
- Maximum flow rate is limited to 10 ml/min .

Fluid type: Viscosity up to 500 cSt

Symbol: H

This setting "Visc500 cSt" is for media with a viscosity of up to approx. 500 cSt. This setting alters the characteristics of the pump as follows:

- The pump suction rate is significantly reduced →
- → The maximum flow rate is limited to 4 ml/min

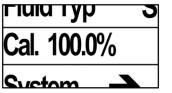


Fig. 41: Calibration

riula ry	p 2
Target val.	20.0 ml/min
Actual val.	20.0 ml/min
Suctor	

Fig. 42: Inputting the calibration actual value

7.12 To calibrate the pump

The pump has already been precisely adjusted in the factory. However, the device will operate even more precisely if it is calibrated. Calibration eliminates any variances caused by supply lines, back-pressure and viscosity.

Calibration can be used to adjust the pump flow rate by $\pm 20\%$.

- 1. Set up system
- 2. Take an accurate measurement of the flow rate or metering volume
- 3. Use the control knob to make the required menu selection [Cal. 100.0%]
- 4. Input the measured actual value with the control knob and confirm (see Fig. 42)
- 5. Check the calibration with a repeat measurement.

NOTE

The calibration function uses the specified flow setting as the target value. This means the pump can always be calibrated for the current operating point.

NOTE

If the target flow rate still cannot be achieved after several calibration attempts, see "Troubleshooting" (Chapter 12).

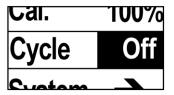


Fig. 43: Cycle display

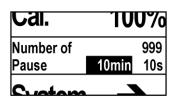


Fig. 44: Setting menu



100.070

Fig. 45: Main display

Udi.

Back

Fnalish

stem

Fig. 46: System menu

Fig. 47: Back to main menu

7.13 Cycle metering mode

The "Cycle" settings in the main menu allow several volumes to be metered with a specific pause time.

- 1. Use control knob to set desired volume
 - Set metering volume as volume in [ml].
 - Set metering volume as flow rate and time in [ml/min].
- 2. Select "Cycle" menu (see Fig. 43)
- 3. Press control knob
 - ➔ The display in Fig. 44 appears
- 4. Turn control knob
 - → Specify number of cycles (Off, 2...999, unlimited)
- 5. Press control knob
- 6. Specify pause time
 - → The following intervals are possible: 1s...99min 59s
- 7. Press control knob

→

Main display appears (see Fig. 45)

NOTE

If the pump is disconnected from the mains or switched off, cycle operation is deactivated for safety reasons.

NOTE

The "PAUSE" button can be used to temporarily interrupt the process. The process can be restarted by pressing the "PAUSE" button again.

7.14 Setting the system

Select "System" to call up a second menu allowing the other pump settings to be entered.

7.15 Back to main menu

Click on "Back" to return to the main menu for entering pump default settings.

Other options for returning to the main display are as follows:

- Press the "STOP" button: this ends the input operation, and the display switches back to the main display
 - ➔ The pump will stop when this operation is carried out
- Press the "START" button: this ends the input operation, and the display switches back to the main display

→ The pump starts when this operation is carried out

- > Wait for 10 seconds without entering any inputs
 - ➔ The pump ends the input operation and switches back to the main display



Fig. 48: Language selection



7.16 Select language setting

The language setting operation selects the required language for the operating menu.

7.17 Activate autostart

Automatic start-up when autostart function is activated		
The pump starts up by itself and without warning.		
Do not switch on the autostart function until the system has been tested and is ready for operation		
Clearly identify the fact that the autostart option for the		

- the pump has been activated
- Before start-up, check that hoses and equipment are \triangleright leak-tight and correctly connected
- Do not operate the pump with hazardous media \triangleright

Autostart and external drive

Οαι	Alam
Auto start	Off
Contract	10

Fig. 49: Autostart



Fig. 50: Autostart symbol in display



Fig. 51: Contrast

If you want the pump to respond automatically to external signals as soon as the power supply is switched on, you can activate the autostart option.

The "Autostart" setting defines the process for switching on the pump.

Input	Meaning
Off	Pump does not start by itself
	Pump starts automatically when the power supply to the pump is switched on.

Tab. 20: Autostart

When the "Autostart" option is on, this is indicated in the main display with the "AS" symbol (see Fig. Fig. 50).

7.18 Set contrast

Display contrast adjustment. Adjust the contrast if the display is difficult to read.



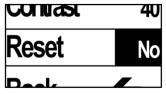


Fig. 52: Reset

7.19 Resetting the pump (Reset)

The "Reset" function is used to reset the pump to its factory settings.

Input	Meaning
No	No reset takes place
Yes	Pump is reset to its factory settings
Code	No function

Tab. 21: Reset

The reset function will restore the following settings:

Value	Setting = factory setting
Flow rate [ml/min]	10.0
Unit	ml/min
Time counter	Offmin sec
Fluid type	Standard
Calibration	100%
Language	English
Analog input	Off
Digital input 1	Off
Digital input 2	Off
Digital output	Alarm
Autostart	Off
Contrast	40

Tab. 22: Reset factory settings

7.20 Shutting down the pump in an emergency

In order to shut the pump down in an emergency, pull out the mains plug. Or shut the pump down using the "Stop" button

RC version (remote control) 8

WARNING

Danger of automatic start-up

The pump starts up by itself and without warning.

- > Do not switch on the autostart function until the system has been tested and is ready for operation
- Clearly identify the fact that the autostart option for the pump has been activated
- Before start-up, check that hoses and equipment are leak-tight and working properly
- Do not operate the pump with hazardous media \geq

These remote control functions are only available in the RC version.

RC cable

- 1. Connect cable for remote control (RC cable) to a suitable signal source.
 - 1. Remove protective caps from RC connector plug.
 - 2. Connect RC cable to pump.
- Analog input 1
- Brown
- 2 **Digital input 1**
 - White
- 3 **Digital input 2**
 - Blue
- **Open collector output** 4
- Black
- Ground 5
 - Grey

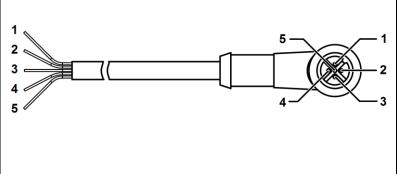
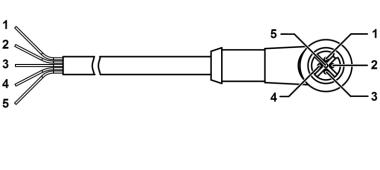


Fig. 53: KNF Flodos RC cable pin assignment

Pin	Wire colour	Description	Function	max. value
1	brown	Analog input	0.15% to 100% up to flow rate, depending on settings	10 V DC
2	white	Digital input 1	Start / stop	24 V DC
3	blue	Digital input 2	Reset / Prime / Foot Switch	24 V DC
4	black	Digital output	Various output signals	open collector max. 10 mA max. 24 V DC
5	grey	Ground		

Tab. 23: RC cable pins



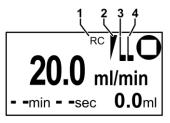


Fig. 54: RC main display

Activate analog input

When external drive is activated, this is indicated in the main display by the "RC" symbol (1) (see Fig. Fig. 54).

According to the inputs activated, the display also shows symbols for analog input (2), digital input 1 (3) and digital input 2 (4).

8.1 Remote control analog input

The analog input can be used to externally specify a pump flow rate (see Tab. 26).

The analog input can only be activated in transfer operation [ml/min].

NOTE

When analog input is active:

- the pump can be started only if a valid analog signal is present
- the time counter is deactivated
- manual inputting of flow metering volume is disabled

A_in Off

Fig. 55: Analog input

Control signals

Input	Meaning	Signal
Off	Analog input not active	
0-5 V	Voltage control [V]	0.05 - 5
0-10 V	Voltage control [V]	0.1 - 10
4-20 mA ⁷	Current control [mA]	4.16 - 20
0-20 mA	Current control [mA]	0.2 - 20

Tab. 24: Analog input options

On/off thresholds

In order to obtain clear switching criteria, a switching hysteresis is implemented.

Input	On threshold	Off threshold
0-5 V	0.05 V	0.04 V
0 – 10 V	0.1 V	0.09 V
4 – 20 mA	4.16 mA	4.08 mA
0 – 20 mA	0.2 mA	0.18 mA

Tab. 25: On/off thresholds

⁷ If the input is less than 2 mA, the pump will go into alarm mode, and the alarm is automatically reset as soon as the input exceeds 2 mA.


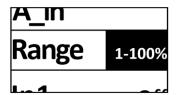


Fig. 56: Flow rate range selection

Flow rate range selection for analog input

There are three flow rate ranges for different set points (see Tab. 26).

•	100%:	1 – 100%	0.2– 20 ml/min
•	30%:	0.3 – 30%	0.06 – 6 ml/min
•	15%:	0.15-15%	0.03 – 3 ml/min

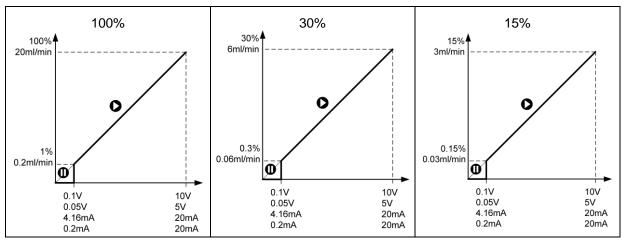

START / STOP analog signal

Tab. 26: Analog Input

Starting and stopping the pump with an analog signal

- 1. Set analog input to correct signal type
- 2. Press the "START" button.
 - → Pump switches to "Pause" condition.
- 3. Apply analog signal.
 - → The pump now starts operating, and switches to "Run" condition.
- 4. Reduce analog signal to below cut-off threshold.
 - → Pump stops and switches to "Pause" condition.

NOTE After the pump has been stopped, there is a pause time of 0.2 seconds before it can be restarted.

Stopping the pump with the "STOP" button is called a "Master Master Stop Stop". The pump can then only be restarted manually (see Section 7.6).

NOTE

The pump is supplied with Master Stop. For initial start-up with external drive, it is therefore essential to press "START" once to cancel Master Stop.

For an external restart after a power cut, use autostart (see Section 7.17) or Reset on digital input 2 (see Section 8.3).

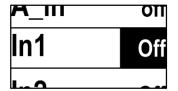


Fig. 57: Digital input 1

8.2 Digital input 1, Start / Stop

The pump can be started and stopped externally via digital input 1.

- If digital input 1 is active, the pump can only be started via the digital input.
- If analog actuation is also used, a valid analog signal must be present.

Control signals

Input	Meaning	Signal
Off	Digital input 1 not active	
Level	Start / Stop according to signal level	< 0.8V = On > 2.0V = Off
Pulse	Start / Stop on signal pulse	< 0.8V = On > 2.0V = Off

Tab. 27: Digital input 1 options

NOTE

To ensure that the digital input operates correctly:

- after the pump has been stopped, there must be a pause time of 0.2 seconds
- > select a pulse duration of more than 0.1 seconds
- use debounced switches

START / STOP in "Pump ml/min without timeout" condition

Start / Stop level

- Set digital input 1 "In_1" to Level.
 Press the "START" button.
 - ➔ Pump switches to "Pause" condition.
- 3. Apply signal 1.
 - → The pump will start to operate if the level is below 0.8 V, and stop if it is above 2.0 V (see Fehler! Verweis-quelle konnte nicht gefunden werden.).

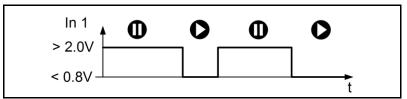


Fig. 58: Start / Stop level actuation

- Start / Stop pulse
- Set digital input 1 "In_1" to Pulse
 Press the "START" button.
 - ➔ Pump switches to "Pause" condition.
- 3. Apply signal 1.
 - ➔ The pump will start to operate when a pulse of less than 0.8V is generated, and stop on the next pulse of more than 2.0V.

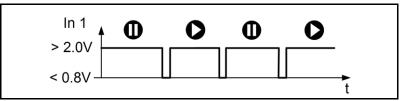


Fig. 59: Start / Stop pulse actuation

Start / Stop level with timeout

Start / Stop level with timeout

- 1. Set digital input 1 "In_1" to Level.
- 2. Press the "START" button.
 - ➔ Pump switches to "Pause" condition.
- 3. Apply signal 1.
 - ➔ The pump will start to operate when the level is connected to ground, and stop when the level rises above 2.0 V.


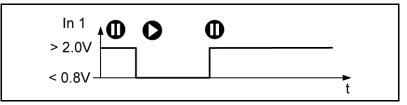


Fig. 60: Start ml/min with time counter level control

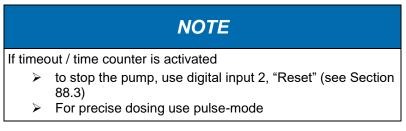

Start / Stop pulse with timeout

- 1. Set digital input 1 "In_1" to Pulse
- 2. Press the "START" button

Pump switches to "Pause" condition.

3. Apply signal 1

→

→ The pump will start to operate when a pulse to ground is generated, and stop when the set time has elapsed.

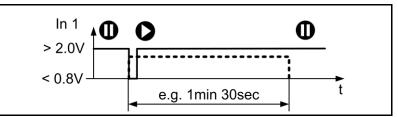


Fig. 61: Start ml/min with time counter pulse control

8.3 Digital input 2 Reset / Prime / Foot Switch

Digital input 2 can be used for an externally activated error "Reset" or fluid system "Priming" operation.

The level and pulse settings are the same as for digital input 1.

Digital input 2 is mainly used with a foot switch (see Section 13.5).

Control signals

Input	Meaning	Signal
Off	Digital input 2 not active	
Reset	Reset of pump errors and cancel metering with time counter	< 0.8V = Reset
Priming	Rapid priming of pump head and hoses (see Section 7.12)	< 0.8V = Prime
Combined	Reset of pump errors and cancel metering with time counter, then prime	< 0.8V reset for falling signal edge Prime after 1 second < 0.8V
Level	Start / Stop according to signal level	< 0.8V = On > 2.0V = Off
Pulse	Start / Stop on signal pulse	< 0.8V = On > 2.0V = Off

Tab. 28: Digital input 2 options

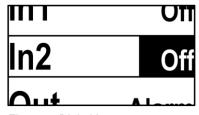


Fig. 62: Digital input 2

Reset Digital input 2 Reset

The "Reset" setting activates the following functions on application of a signal to digital input 2:

- Interrupt a metering process
- Reset pump errors
- Cancellation of an active Master Stop, pump is now in "Pause" operation condition, and responds to external control signals.

Priming Digital input 2 Prime

The "Prime" setting activates the following functions on application of a signal to digital input 2:

- Interrupt a metering process
- Reset pump errors
- Pump runs in high-speed "Prime" mode for the duration of the signal
- Cancellation of an active Master Stop, pump is now in "Pause" operation condition, and responds to external control signals.

Combined **Digital input 2 Combined**

The "Combined" setting activates the following functions on application of a signal to digital input 2:

- Interrupt a metering process
- Reset of pump errors on signal edge
- If the signal is applied for longer than 1 second, the pump will operate in "Prime" at high speed for the duration of the signal
- Cancellation of an active Master Stop, pump is now in "Pause" operation condition, and responds to external control signals.

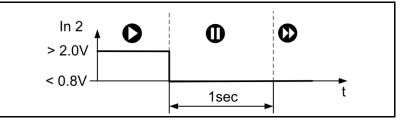


Fig. 63: Combined function sequence

Start / Stop level

START / STOP with foot switch using level

- 1. Set digital input 2 "In_2" to Level.
- 2. Connect foot switch to pump.
- 3. Press "START" the pump goes into "Pause" condition.
 - The pump will keep operating as long as the foot switch is pressed.

NOTE

If time counter is switched on, pump will perform a complete metering cycle.

Start / Stop pulse

top pulse START / STOP with foot switch using pulse

- 1. Set digital input 2 "In_2" to Level.
- 2. Connect foot switch to pump.
- 3. Press "START" the pump goes into "Pause" condition.
 - → Each time the foot switch is pressed the pump is started or stopped.

NOTE

If timeout / time counter is switched on, pump will perform a complete metering cycle.

8.4 Digital output

The digital output can be used to feed back a number of messages to the control system.

The output is an open collector output (see Section 3.8).

Output signals

Set output function to:	Meaning	Signal
Alarm	Pump error signal	Level
Motor	Signal when pump motor running	Level
Vol.End	Signal when volume metering ends	Level
Mot.Pul.	10 pulses per revolution of the pump motor	Pulse
Vol.Pul.	One pulse per 20 μ l of fluid transferred	Pulse

Tab. 29: Digital output options



9 RC Plus version (serial communication)

A WARNING

Danger of automatic start-up

The pump starts up by itself and without warning.

- Do not transmit a start command until the system has been tested and is ready for operation
- Mark remote-controlled pumps
- Before start-up, check that hoses and equipment are leak-tight and working properly
- > Do not operate the pump with hazardous media

The functions for remote control via the RS 232 interface are only available with the RC Plus version.

RC cable

- 1. Connect cable for remote control (RC cable) to a suitable signal source.
 - 2. Remove protective caps from RC connector plug.
 - 3. Connect the RC cable to the pump.
 - 4. The commands and protocol are available on <u>www.knf.com/downloads</u> (search: "communication protocol")

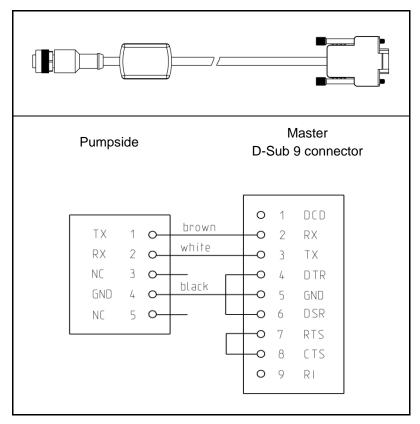


Fig. 65: KNF Flodos RC cable pin assignment



10 Cleaning and maintenance

10.1 Safety

Danger of injuries caused by hazardous materials

Poisoning and caustic burns caused by contact with aggressive, inflammable or radioactive media

- > Flush the pump thoroughly (see Chapter 10.2)
- Decontaminate the pump if necessary
- Observe the safety data sheets for the media to be transferred
- > Wear suitable protective equipment

A WARNING

Gefahr vor Wärmeentwicklung durch chemische Reaktionen mit der Spülflüssigkeit

Starke Wärmeentwicklung schädigt die Membrane, O-Ringe und das Ventilsystem, dies führt zu Leckagen

- > Chemische Reaktionen in der Pumpe vermeiden
- Sicherheitsdatenblätter der geförderten Flüssigkeit beachten
- Nachfolgende Ratschläge beachten

The following requirements apply to all the activities described in this chapter:

- System depressurised
- Specialist personnel
- The type of protective equipment required will depend on the medium to be transferred.
- Observe the safety data sheets for the media to be transferred.

Prior requirements Personnel Personal protective equipment

KNF Flodos | BA_SIMDOS02_EN_09_166267 | Translation of original operating instructions, English

10.2 Cleaning the pump

In order to maintain the service life of the pump, after each use and before any long periods of inactivity the pump must be flushed through with neutral media.

IMPORTANT:

To ensure satisfactory start-up when the unit is again required, it is important to ensure that the pump is free of any crystallising, adhesive or curing media.

10.2.1 Metering medium: acids

- Flush pump with a suitable neutralising liquid in a circular flow for 10 minutes
- > Then flush with distilled water for 5 minutes

10.2.2 Metering medium: bases

- Flush pump with a suitable neutralising liquid in a circular flow for 10 minutes
- > Then flush with distilled water for 5 minutes

10.2.3 Metering medium: organic solvents

 Flush pump with isopropyl alcohol (C3H8O) in a circular flow for 10 minutes

NOTE

Never flush non-polar solvents immediately with water

10.2.4 Metering medium: biological solution

- Flush pump with 10% hydrogen peroxide in a circular flow for 10 minutes
- > Then flush with distilled water for 10 minutes




The procedure described in Section 10.2.4 is not a sterilisation

Interval	Assembly	Action	Personnel
When changing the transfer medium	Complete pump	5. Flushing the pump	Specialist per- sonnel
Daily	Complete pump	 Visually inspect the pump for pollution/contamination. If pollution/contamination is found, clean the pump. Check the discharge outlet on the intermediate flange for leaks 	Trained person- nel
	Complete pump	 9. Visually inspect the pump for external damage. 10. If damage is found, take corrective action (see Chapter "Troubleshooting") 	Trained person- nel
	Complete pump	11. Visually inspect the pump for leaks12. If leaks are found the pump must under no circumstances be used. (see "Troubleshooting")	Specialist per- sonnel
	Complete pump	 13. Check pump capacity 14. If pump capacity falls, clean the pump head and replace elastomer parts where necessary. 	Specialist per- sonnel
Weekly	Filter (accessory)	15. Check filter for contamination.16. If contamination is found, clean or replace the filter	Specialist per- sonnel
1,000 hours of operation	Pump diaphragm, valves and seal	17. Replace pump diaphragm, valves and seal	Specialist per- sonnel
as required	Pump head	18. Replace pump head	Specialist per- sonnel

10.3 Maintenance plan

Tab. 30: Servicing schedule

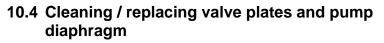

	Poisoning escaping h nance/serv > Fo > En (pa > Ch	om hazardous materials and caustic burns or undesirable reactions caused by nazardous materials as a result of incorrect mainte- vicing or the transferring of incompatible media. Illow the maintenance instructions usure cleanliness during maintenance articles on seal faces cause leaks) neck that pump and system are leak-tight after mainte- nce
	If they com cause pois > Flu pu > We	poisoning and caustic burns ne into contact with skin and eyes, aggressive media soning and caustic burns. Ush pump as described in Section 10.2, Cleaning the mp. ear protective equipment, e.g. protective gloves, safety asses.
Prior requirements	■ Th	imp switched off and disconnected from mains the pump must be free of any hazardous substances pases must be disconnected from the pump head
Personnel Personal protective equipment	SpTh	becialist personnel the type of protective equipment required will depend on the medium to be transferred.
	 Ot 	e medium to be transferred. oserve the safety data sheets for the media to be trans- red.
Tools and materials	Qty	Tools/materials

als	Qty	Tools/materials
	1	Phillips screwdriver no. 1
	1	Spare parts kit (see Section 13.1)
	1	Torque screwdriver that can be set to 0.55 Nm.

Tab. 31: Tools/materials

Information on procedure

Always replace valve disks, valve seats and seals at the same time, to maintain satisfactory flow rate performance.



KNF Flodos | BA_SIMDOS02_EN_09_166267 | Translation of original operating instructions, English

10.4.1 Cleaning the pump head

The pump head should only be cleaned if:

- > the pump is no longer sucking properly,
- > the flow rate is too low,
- > reverse suction can be seen, or
- the pump head cannot be cleaned satisfactorily by flushing.

NOTE

Ensure cleanliness when replacing valves, as pollution/contamination may prevent the valve from working properly.


- 1. Disassemble the pump head as described in Section 10.4.2
- 2. Rinse connecting plate (3) and intermediate plate (8) and wipe with a soft cloth. Blow dry with compressed air
- 3. Remove valve seats (5) carefully. The sealing washers (4) are not present in the FT head, see Fig. 66
- 4. Remove O-Rings (6) and valve disk (7).

NOTE

Do not damage the sealing rings.

- Do not use any abrasive agents.
- > Do not clean head parts with hard materials.
- Check head parts, especially sealing edges and seal faces, for damage
- Replace any damaged parts
- 5. Wash all parts with flushing agent and blow dry with compressed air
- 6. Wipe the pump diaphragm with a soft damp cloth
- 7. Assemble the pump head as described in Section 10.4.2
- 8. Check that the pump is leak-tight (see Section 10.5)
 - 1 2 3 4 5 6 7 8

Fig. 66: Head, KT, TT

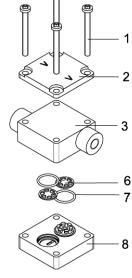

Head, FT


- 1 Head screws
- 2 Head plate
- 3 Connecting plate
- 4 Sealing washer
- 5 Valve seat
- 6 O-ring
- 7 Valve disk
- 8 Intermediate plate

10.4.2 Replacing the valve




> marking if it deviates from the standard

Ensure cleanliness when replacing valves, as pollution/contamination may prevent the valve from working properly.

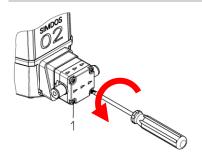


1. Undo head screws (1)


- Completely remove head (10)
 Remove head plate (2) and head screws (1)
 - 4. Check diaphragms (20) for damage and replace where necessary

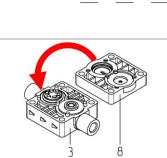

5. Fold the connecting plate (3) and the intermediate plate (8) apart

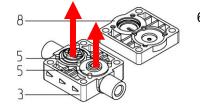


6. Remove valve seats (5) carefully

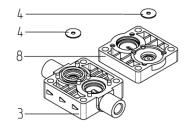

Replace sealing washers (4)
 The FT head does not contain a sealing washer or valve seats

Caution!

Sealing washer must lie flat in the valve seat, otherwise pump will not be leak-tight

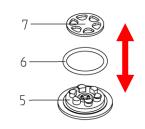


- 8. Remove valve disks (7) and O-rings (6)
- 9. Fit O-rings (6) flat on the base, taking care not to twist them
- 10. Fit valve disks (7), which must lie flat on the valve

Caution! Observe material coding (below)

11. Check valve material before using. For further information, see Section 4.3




DuPont™Kalrez®

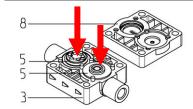


12. Place pre-assembled valve seats (5) in connecting plate (3)




CHEMRAZ®

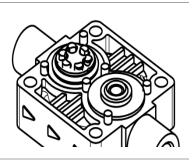

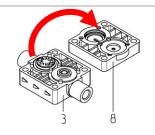

13. Fit the connecting plate (2) and the intermediate plate (1) together

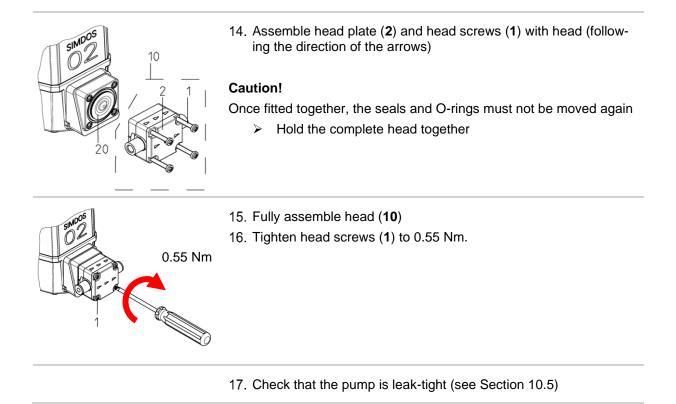

10.4.3 Replacing the pump diaphragm

- Prior requirements
 Pump must be free of any hazardous substances
 - Hoses must be disconnected from the pump head
 - Pump is connected to the mains
 - specialist personnel trained by KNF
 - The type of protective equipment required will depend on the medium to be transferred.
 - Observe the safety data sheets for the media to be transferred.

Ensure cleanliness when replacing diaphragms, as pollution/contamination may prevent the pump from working properly.

- 1 Support / Connecting rod
- 2 Insert ring
- 3 Pump diaphragm

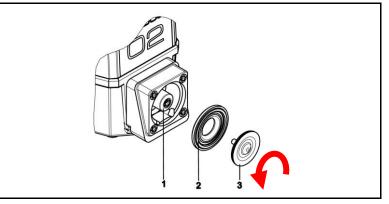


Fig. 67: Replacing the pump diaphragm

Removal

- 1. Activate the diaphragm maintenance function in the pump system menu
- 2. Disassemble pump head as described in Section 10.4.2
- 3. Lift pump membrane (**3**) at opposite side edges, grip and screw out by turning anticlockwise (see arrow in Fig. 67)
- 4. Inspect all parts for pollution/contamination and damage, cleaning them if necessary

Installation

- 5. Activate diaphragm maintenance in the pump system menu (if no longer active)
- 6. Fit insert ring (2).

→

- 7. Screw pump membrane (3) clockwise onto connecting rod (1) and hand-tighten.
- 8. Press the "STOP" button,
- 9. Activate "Prime" in main menu for 2s
 - Diaphragm moves into housing
- 10. Assemble the pump head as described in Section 10.4.2.
- 11. Check that the pump is leak-tight (see Section 10.5)

NOTE

If the pump diaphragm is not fitted correctly, the pump may be mechanically obstructed and become damaged as a result.

- Screw the pump membrane in properly as far as it will go
- Do not damage the surface of the diaphragm

Personnel Personal protective equipment

10.5 Checking that pump is leak-tight

Danger of bursting of the fluid system as a consequence of overpressure.

The pump builds up pressure. In a closed system the max. permitted operating pressure may be exceeded, thus causing damage to the pump and/or the system.

- In the case of parts that are in contact with the fluid, only select those that are designed for the pump's operating pressure
- Do not transfer against shut-off devices or closed systems, or
- install pressure relief/safety valves and set them to 6 bar (in the KNF product range)

In order to ensure correct assembly and hence reliability, it is essential to perform a leak test.

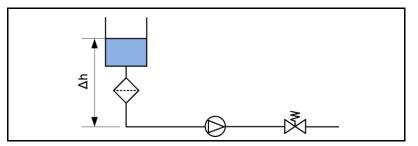


Fig. 68: Pressure test setup

- 1. Set up the system as shown in Fig. 68.
- 2. Use distilled water as test medium.
- 3. Set pressure control valve to working pressure.
- 4. Make sure that the fitting is leak-tight (see Section 6.2.3).
- 5. Operate pump carefully for 5 minutes.
- 6. Check pump for leaks.

If no traces of medium are visible on the head, the pump can be used in accordance with its specification (see Chapter 3).

Caution!

If it is found that liquid is escaping, under no circumstances should the pump continue to be used.

- 1. Check that connections are leak-tight.
- 2. Check tightening torque of head screws.
- 3. Check that elastomer parts are clean.
- 4. Check leak-tightness again.

If these actions have no effect, contact your local KNF dealer and do not under any circumstances continue to use the pump.



11 Shutdown

11.1 Safety

A DANGER

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
Danger of injuries caused by hazardous materials

Poisoning and caustic burns caused by contact with aggressive, inflammable or radioactive media

- Flush the pump thoroughly (see Chapter 10.2)
- Decontaminate the pump if necessary
- Observe the safety data sheets for the media to be transferred
- > Wear suitable protective equipment

The following requirements apply to all the activities described in this chapter:

- System depressurised
- Specialist personnel
- The type of protective equipment required will depend on the medium to be transferred.
- Observe the safety data sheets for the media to be transferred.
- 1. Flush the whole system including the pump with a neutral liquid
- 2. Pump the pump empty (see Section 10.2).
- 3. Press the STOP button to end the transferring operation
- 4. Unplug the pump from the power supply
- 5. Dispose of pump properly



Prior requirements

Personnel

Personal protective equipment



12 Troubleshooting

12.1 Safety

none

•

Disconnect the mains plug from the power supply before working on the pump head.

Prior requirements

Personnel

Pump has been thoroughly flushed/decontaminated Specialist personnel

Personal protective equipment

12.2 Displaying of error messages

Display	Description	Fault remedy
Error 1 Motor	Control deviation too high, motor is over- loaded	Switch pump on / off
Error 2 Temperature	Motor overheating	 Allow pump to cool Reduce ambient temperature
Error 3 Supply	Supply voltage is less than 21 V	 Supply with 24 V and sufficient power
Error 4 Encoder	Position measuring malfunction	Switch pump on / off
Error 5 4 – 20 mA	Analog set point setting less than 2 mA	Check control signalCheck cable
Error 6 Flash	Error in memory	Switch pump on / off
Error 7 Overpressure	System pressure exceeds 7 bar	 Check pump for closed valves and blocked fil- ters

If the fault cannot be rectified, please contact your local KNF dealer (see last page for address)

12.3 Rectifying	transfer problems	
Symptom	Possible cause	Fault remedy
Pump cannot be switched on	Pump not connected to the mains.	 Connect pump to the mains. Check mains voltage
Pump does not work	Connections or hoses are blocked.	 Check hoses and connections. Remove blockage.
Pump does not work	External valve is closed or filter is clogged or too small.	1. Check external valves and filters.
Flow rate too low	Particles in the pump head	 Clean valves (see Maintenance) Clean pump head (see Maintenance)
Insufficient vacuum		 Use prefilter
Flow rate too low	Materials chemically dam- aged by pumped media.	 Select a type of material that is resistant and suitable (see chemical resistance list).
		2. Replace valves (see Section 10.4)
Flow rate too low	Worn diaphragm or valve plates/seals.	 Replace diaphragm and valve plates/seals (see Section 10.4).
Pump not drawing any vacuum	Pump can only draw vacu- um against atmosphere.	 Change pressure conditions on pres- sure side.
Unstable flow rate	Cross-section of hydraulic hoses or connectors too	 Disconnect the pump from the system and determine output values.
	narrow or restricted.	 Remove restriction (e.g. valve) if neces- sary.
		If applicable, use larger-diameter hoses or connectors.
Leakage in pump	Worn diaphragm or valve	1. Tighten head screws to 0.55 Nm
head	plates/seals.	 Replace diaphragm and valve plates/seals (see Section 10.4).
Leakage from dis- charge port.	Leakage at pump diaphragm	 Contact KNF dealer Send pump to KNF dealer for safety check

12.3 Rectifying transfer problems

Tab. 32: Pump does not work

If the fault cannot be rectified, please contact your local KNF dealer (see last page for address)

Returning the pump

- 1. Flush the pump to clear the pump head of any hazardous or aggressive liquids.
- 2. Disconnect the pump.
- 3. Send the pump, with completed decontamination declaration (see Chapter 14), to KNF Customer Service stating the nature of the pumped medium (see Arranging for return).

13 Spare parts and accessories

13.1 Spare parts

Spare part		Order No.
SIMDOS 02 KT Complete Head Kit;	Standard Kalrez® (without diaphragm)	166297
SIMDOS 02 TT Complete Head Kit;	Standard Kalrez® (without diaphragm)	166298
SIMDOS 02 FT Complete Head Kit;	Standard Kalrez® (without diaphragm)	164981
SIMDOS 02 FT Complete Head Kit;	Optional, Chemraz® (without dia-phragm)	164982
SIMDOS 02 ST Complete Head Kit;	Standard Kalrez® (without diaphragm)	166299
SIMDOS 02 ST Complete Head Kit;	Optional, Chemraz® (without diaphragm)	166300
SIMDOS Valve Kit	Standard Kalrez®	164983
SIMDOS Valve Kit	Optional, Chemraz®	164984
SIMDOS 02 Diaphragm Kit	Standard PTFE	165442
SIMDOS 02 Diaphragm Kit	Standard FFKM	170867

13.2 Hoses

Hoses		Order No.
Connection set, AD1/8, UNF 1/4", L=1m;	(2 Ferrules and 2 Fittings)	069790
Hose, FEP, UNF 1/4", screw connection L=1m	(1 Fitting)	166329
Hose, FEP, UNF 1/4", screw connection L=2m	(1 Fitting)	166330
Hose, FEP 1/16", Gripper, UNF 1/4", L=1m	(1 Fitting)	166335

13.3 Pressure control valve

Pressure control valve - kit	Order No.
SIMDOS 02 Accessory Set JFDV 30 KT	166286
SIMDOS 02 Accessory Set JFDV 30 TT	166287
SIMDOS 02 Accessory Set JFDV 30 FT	166288

13.4 Filter

Included Filter	Head material	Mesh size	Material*	Order No.
Filter FS 60 T	ТТ	70µm	PVDF	165210
Filter FS 60 X	KT, FT, ST	35µm	PEEK	323625 (1 pc)
Filter FS 60 X	KT, FT, ST	35µm	PEEK	323626 (10 pcs)

*material for housing and filter mesh


13.5 Foot switch

Foot switch	Order No.
Pulse foot switch	155872

13.6 Fastening material

Fasteners	Order No.
Tripod holder	160474
Mounting plate	160473

14 Decontamination declaration

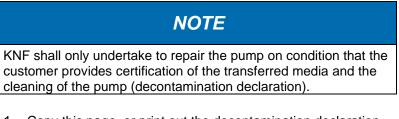

- 1. Copy this page, or print out the decontamination declaration from our website <u>http://www.knf.com/downloads</u>.
- 2. Enter the pump model, the Serial No. and the transferred media in the form below and send the signed form together with the flushed and cleaned pump to KNF Customer Service.

Customer decontamination declaration for repair order

We confirm that the pump below has been used to pump the following media, and that the pump has been flushed and cleaned.

Pump model	
Serial No.	
Pumped media	

The pump does not contain aggressive, biological, radioactive, poisonous, or other dangerous media.

Company

Date/Signature

KNF Flodos AG Wassermatte 2 6210 Sursee, Switzerland Tel +41 (0)41 925 00 25 Fax +41 (0)41 925 00 35 www.knf.com



www.knf.com