


OEM

N87.9
TRANSLATION OF ORIGINAL OPERATING AND
INSTALLATION INSTRUCTION
ENGLISH

DIAPHRAGM PUMP

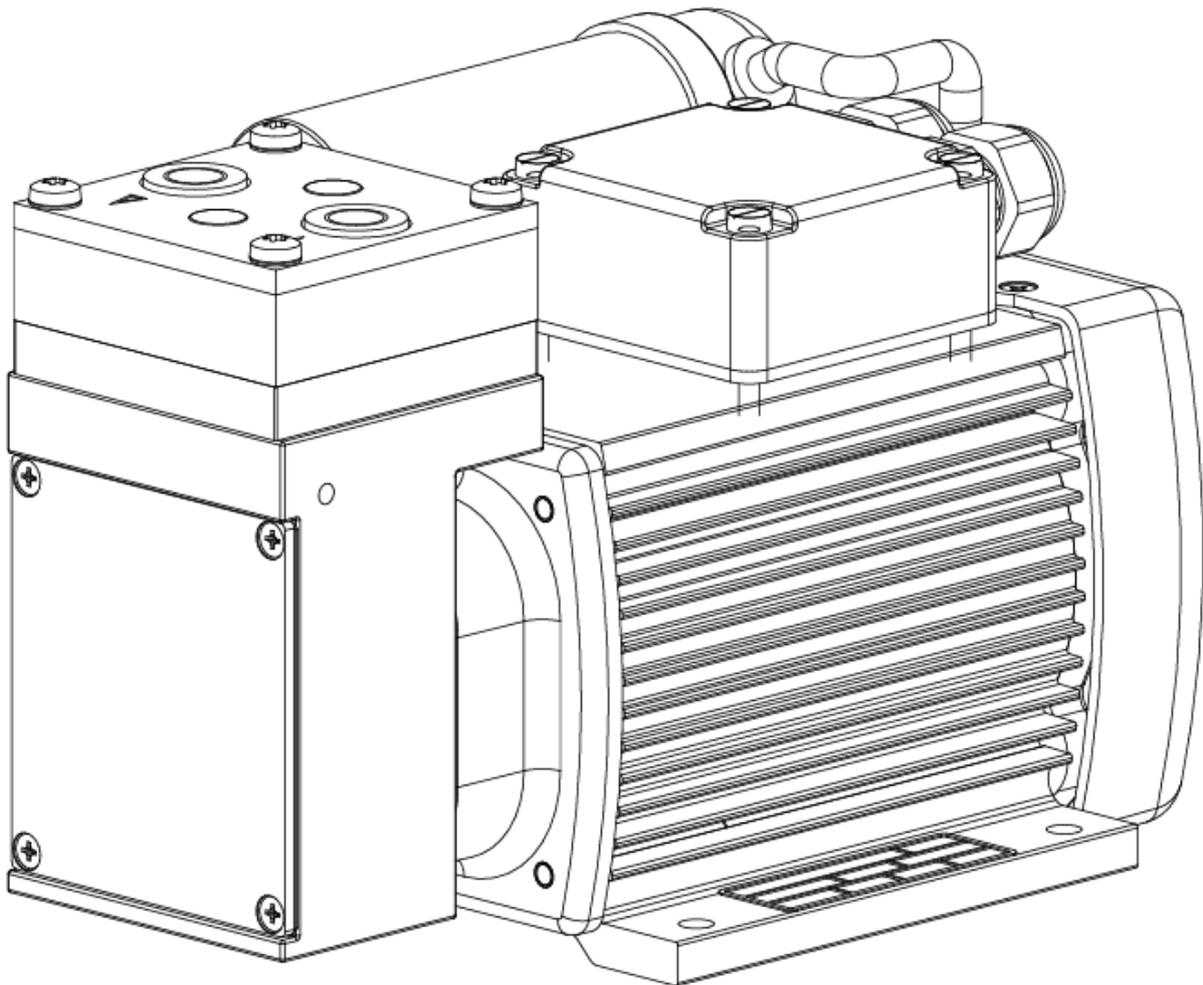

Notice!

Before operating the pump and accessories, read and observe the operating and installation instructions as well as the safety information!

KNF Neuberger GmbH
Alter Weg 3
79112 Freiburg
Germany
Tel. +49 (0)7664/5909-0
Fax. +49 (0)7664/5909-99

www.knf.com

Index

1	About this document	3
2	Use	6
3	Safety	7
4	Technical data	9
5	Design and function	11
6	Transport	13
7	Installation and connection	14
8	Operation	17
9	Servicing	20
10	Troubleshooting	27
11	Spare parts and accessories	29
12	Returns	30

1 About this document

1.1 Using the operating and installation instructions

The operating and installation instructions are part of the pump.

- In the event of uncertainties with regard to the content of the operating and installation instructions, please contact the manufacturer (contact data: see www.knf.com). Please have the type and serial number of the pump ready.
- Read the operating and installation instructions before you commission the pump.
- Give the operating and installation instructions only completely and unchanged to the next owner.
- Keep the operating and installation instructions within reach at all times.

Project pumps For customer-specific project pumps (pump models that begin with "PJ" or "PM"), there may be deviations from the operating and installation instructions.

- For project pumps, also observe the agreed specifications.

1.2 Exclusion of liability

The manufacturer assumes no liability for damages and malfunctions resulting from failure to observe the operating and installation instructions.

The manufacturer assumes no liability for damages and malfunctions resulting from changes or modifications to the device and improper handling.

The manufacturer assumes no liability for damages and malfunctions resulting from impermissible spare parts and accessories.

1.3 Symbols and markings

Warning notice



A notice that warns you of danger is located here.

Possible consequences of a failure to observe the warning notice are specified here. The signal word, e.g., warning, indicates the danger level.

→ Measures for avoiding the danger and its consequences are specified here.

Danger levels

Signal word	Meaning	Consequences if not observed
DANGER	warns of immediate danger	Death or serious injury or serious damage will result.
WARNING	warns of possible danger	Death, serious injury or serious damage is possible.
CAUTION	warns of a possibly dangerous situation	Minor injury or damage is possible.
NOTICE	Warns of possible damage	Damage is possible.

Tab.1 Danger levels

Other notices and symbols

→ An activity to be carried out is specified here (a step).

1. The first step of an activity to be carried out is specified here.
Follow other sequentially numbered steps.

i This symbol indicates important information.

Explanations of pictograms

Pictogram	Meaning
	General warning symbol
	Warning of electrical voltage
	Warning of poisonous substances
	Warning of hand injuries through crushing
	Observe the operating instructions
	General mandatory sign
	Environmentally conscious disposal

Tab.2 Explanations of pictograms

2 Use

2.1 Proper use

The pumps are intended exclusively for transferring gases and vapors.

Responsibility of the owner

Operating parameters and conditions

Only install and operate the pumps in accordance with the operating parameters and conditions described in Chapter 4 *Technical data*.

Only pumps that are fully assembled and in the condition as delivered may be operated.

Make sure that the installation location is dry and that the pump is protected against rain, splash, gushing, and drip water as well as from other contaminants.

Check the tightness of the connections between the pipes of the application and the pump (or the connection of the pump) at regular intervals. Leaky connections carry the risk of releasing dangerous gases and vapors from the pump system.

Requirements on the transferred medium

Before transferring a medium, check whether the medium can be transferred danger-free in the specific application.

Before using a medium, check the compatibility of the media-contacting components (see 4 *Technical data*) with the medium.

Risk of dangerous gas mixtures during pump operation if diaphragm breaks: Depending on the medium being transferred, breakage of the diaphragm can result in a dangerous mixture if the medium mixes with the air in the compressor housing or the surroundings.

Only transfer gases that remain stable under the pressures and temperatures that arise in the pump.

2.2 Improper use

The pumps are not allowed to be operated in explosive atmospheres.

The pumps are not suitable for transferring:

- Dusts
- Liquids
- Aerosols
- Biological and microbiological substances
- Fuel
- Explosives and flammable materials
- Fibers
- Oxidants
- Food

Pumps that can produce both vacuum as well as overpressure may not be used to simultaneously produce vacuum and overpressure.

This function can be made possible on a project basis following consultation with KNF Customer Service.

No overpressure may be applied to the suction side of the pump.

This function can be made possible on a project basis following consultation with KNF Customer Service.

Pumps with capacitor motor are not intended for operation with a frequency converter.

3 Safety

i Observe the safety notices in Chapters *7 Installation and connection* and *8.1 Operation*.

The pumps are produced in accordance with the generally recognized rules of engineering, as well as the occupational health, safety and accident prevention regulations. Nevertheless, dangers can arise during their use that lead to injuries to the user or third parties or to damage to the pump or other property.

Only use the pumps in perfect technical condition, for their intended use, safely and aware of the dangers and in observation of the operating and installation instructions.

The components that are to be connected to the pumps must be designed according to the pneumatic data of the pumps.

When connecting the pumps to the electrical power, observe the corresponding safety rules.

Personnel Make sure that only specially trained and instructed personnel work on the pumps. This applies, in particular, to assembly, connection and service work.

Make sure that the personnel have read and understood the operating and installation instructions, particularly the chapter on safety.

Working in a safety conscious manner Observe the regulations on accident prevention and safety during all work on the pumps and during operation.

Avoid contact with the heads and housing parts, as the pump heats up during operation and may remain hot for some time after operating.

When working on the pump, make certain that the pump is disconnected from mains and without power.

Ensure that no hazards arise from gas flowing when gas connections are open, from the effects of noise or from hot, corrosive, dangerous and environmentally hazardous gases.

Make sure that an EMC-compliant installation of the pump is ensured at all times to prevent the occurrence of dangerous situations.

Handling of hazardous media Upon breakage of the diaphragm and/or leaks, the transferred medium mixes with the air in the surroundings and/or in the pump housing. Make sure that a dangerous situation cannot arise as a result.

When pumping hazardous media, observe the safety regulations for the handling of said media.

Handling of combustible media Note that the pumps are not designed to be explosion-proof.

Make certain that the temperature of the medium is always sufficiently below the ignition temperature of the medium so as to prevent ignition or explosion. This also applies for abnormal operating situations.

Note that the temperature of the medium increases when the pump compresses the medium.

Therefore, make certain that the temperature of the medium also remains sufficiently below the ignition temperature of the medium even when it is compressed to the maximum permissible operating pressure of the pump. The maximum permissible operating pressure of the pump is stated in Chapter *4 Technical data*.

Make certain that the permissible ambient temperature (*4 Technical data*) is not exceeded.

Where applicable, also take into account external energy sources (such as radiation sources) that could additionally heat the medium.

In case of doubt, contact KNF Customer Service.

Environmental protection	Store and dispose of all replacement parts in accordance with the environmental regulations. Observe the respective national and international regulations. This applies in particular to parts that are contaminated with toxic substances.
EU/EC directives/standards	<p>With respect to the Machinery Directive 2006/42/EC, the pumps are partly completed machinery and are, therefore, to be regarded as not ready for use. Partly completed machinery may not be commissioned until it has been determined that the machine into which the partly completed machinery is to be installed complies with the provisions of the Machinery Directive 2006/42/EC. The following fundamental requirements of Annex I of Directive 2006/42/EC (general principles) are applied and observed:</p> <ul style="list-style-type: none">▪ General principles no. 1 <p>As these partly completed machines are built-in devices, the mains connections and equipment for disconnecting and switching off the partly completed machinery as well as overcurrent and overload protection gear must be considered when mounting.</p> <p>Furthermore, protection against contact with moving and hot parts, if present, must be provided during installation.</p> <p> The pumps comply with Directive 2011/65/EU.</p> <p>The following harmonized standards are satisfied:</p> <ul style="list-style-type: none">▪ EN 55014-1/2▪ EN 60204-1▪ EN IEC 63000
Customer service and repairs	<p>The pumps are maintenance-free. However, KNF recommends periodic inspection of the pump for obvious changes in noise or vibration.</p> <p>Only have repairs to the pumps performed by qualified KNF personnel.</p> <p>Housings with electrically live components may only be opened by specialist personnel.</p> <p>Use only genuine spare parts from KNF when performing servicing work.</p>

4 Technical data

4.1 Technical data

Pump materials

Assembly	Material
Pump head	PVDF
Diaphragm	PTFE-coated
Valves	FFPM

Tab.3 Pump materials

Pneumatic values

Parameter	Value
Max. permissible operating pressure [bar rel*]	1.5
Ultimate vacuum [mbar abs.]	140
Flow rate at atm. pressure [l/min]**	7.5

Tab.4 Pneumatic values

*Bar rel related to 1013 hPa

**Liters in standard state (1013 hPa, 20°C)

Pneumatic connections

Pump type	Value
N87TTE	Thread size G 1/8*

Tab.5 *Acc. to ISO 228

Electrical data

Parameter	Value	Value	Value	Value
Voltage [V]	230	115	100	220
Frequency [Hz]	50	60	50/60	60
Power P ₁ [W]	70	70	70	70
Voltage tolerance	± 10%			
Motor protection class	IP54			
Rated current consumption [A]	0.45	0.7	1.0	0.5

Tab.6

Weight

Pump type	Value [kg]
N87TTE	3.1

Tab.7 Weight

Other parameters

Parameter	Value
Permissible ambient temperature [°C]	+ 5 to + 40
Permissible media temperature [°C]	+ 5 to + 40
Dimensions	See Fig. 3 Chapter 7.1 <i>Installing the pump</i>
Highest permissible relative air humidity of the environment	80% for temperatures up to 31°C, decreasing linearly to 50% at 40°C.
Maximum installation altitude [m above sea level]	2000
Gas tightness* of the pump head (leak rate)	$\leq 6 \times 10^{-3}$ mbar l/s

Tab.8 Other parameters

* The gas tightness of the pump head is no longer ensured after the pump head is opened or after replacing diaphragms and valve plates. A leak test can be used to determine whether the original gas tightness is again achieved.

5 Design and function

Design

- 1 Pump outlet
- 2 Pump inlet
- 3 Pump head
- 4 Electrical terminal box
- 5 Motor

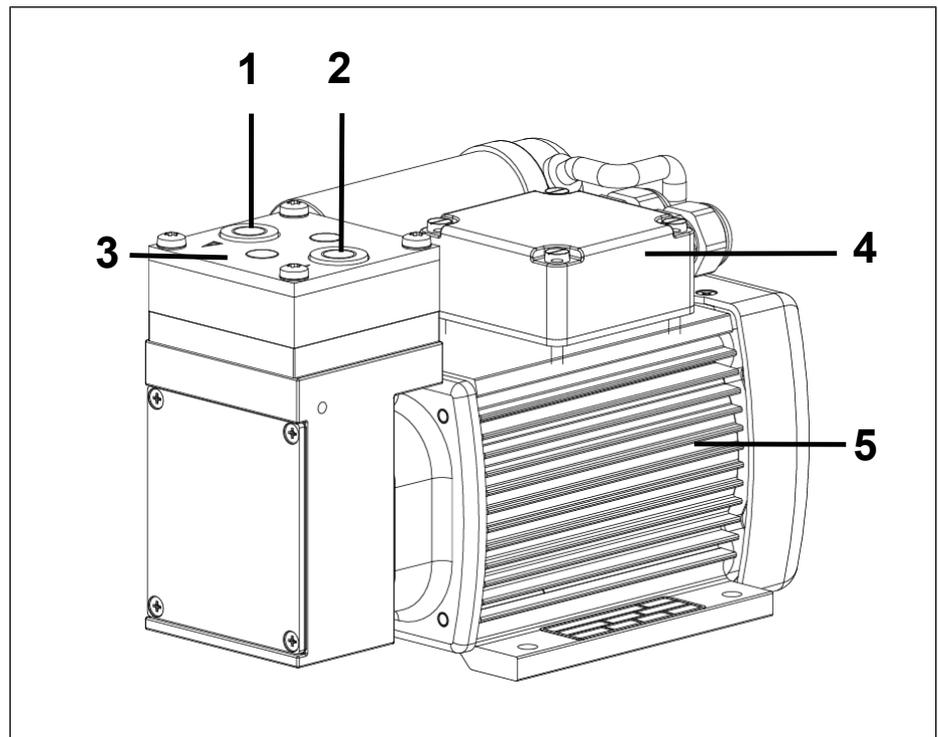


Fig.1 Design N87

Function of a diaphragm pump

- 1 Outlet valve
- 2 Inlet valve
- 3 Transfer chamber
- 4 Diaphragm
- 5 Eccentric
- 6 Connecting rod

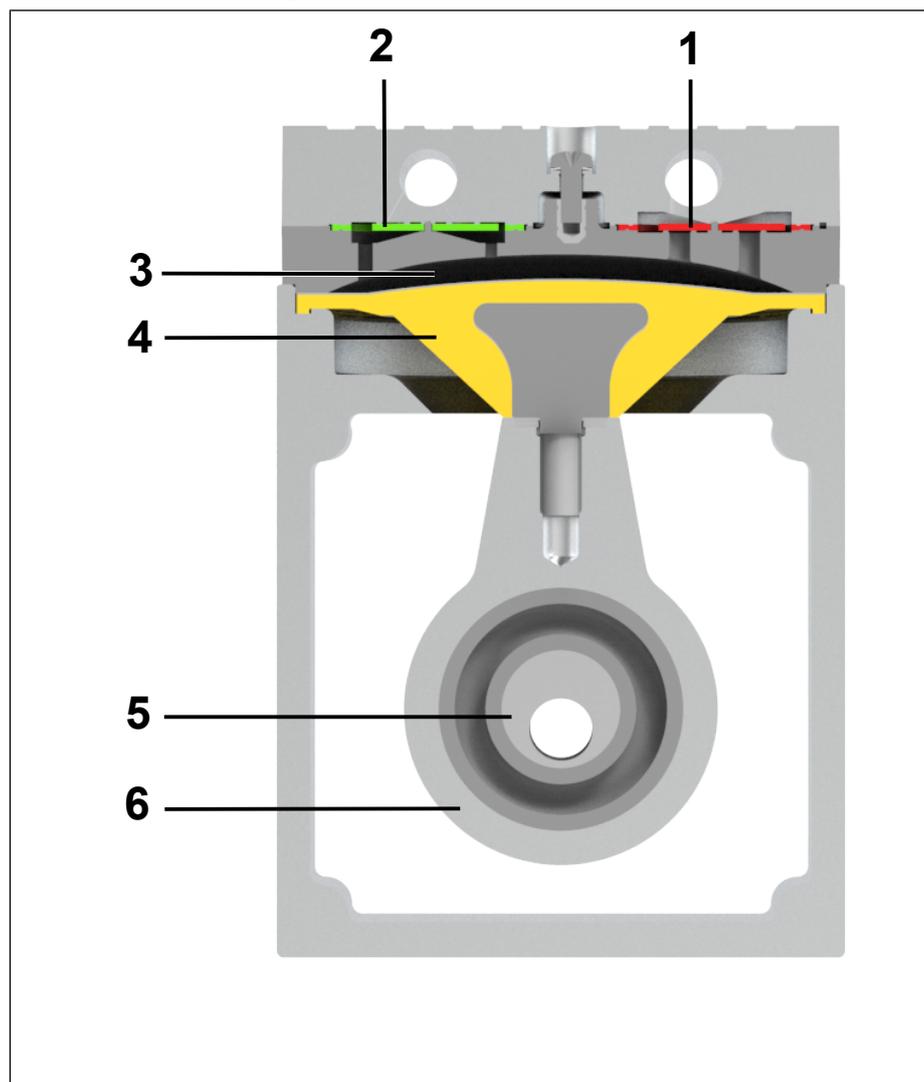


Fig.2 Function of a diaphragm pump

Diaphragm pumps transfer, compress (depending on the version) and evacuate gases and vapors.

The elastic diaphragm (4) is moved up and down by the eccentric (5) and the connecting rod (6). In the downwards stroke, it aspirates the gas to be transferred via the inlet valve (2). In the upwards stroke, the diaphragm presses the medium out of the pump head via the outlet valve (1). The pumping chamber (3) is separated from the pump drive by the diaphragm.

6 Transport

General



Personal injury and/or property damage due to incorrect or improper transport of the pump

In the event of incorrect or improper transport, the pump can fall down, be damaged or injure persons.

- Use suitable auxiliary means if necessary (carrying strap, lifting gear, etc.).
- Where appropriate, wear suitable personal protective equipment (e.g., safety shoes, safety gloves).



Risk of injury from sharp edges on the packaging

There is a risk of injury from cutting on the sharp edges when grabbing corners or when opening the packaging.

- Where appropriate, wear suitable personal protective equipment (e.g., safety shoes, safety gloves).

- Transport the pump in the original packaging to the installation location.
- Store the original packaging of the pump (e.g., for later storage).
- Inspect the pump for transport damage after receiving it.
- Document any transport damage in writing.
- Remove any transport safeguards on the pump prior to commissioning.

Parameter

Parameter	Value
Storage temperature[°C]	+ 5 to + 40
Transport temperature [°C]	- 10 to + 60
Permissible humidity (non-condensing) [%]	30 to 85

Tab.9 Transport parameters



Prior to commissioning, make sure that the pump has reached the ambient temperature (4 *Technical data*).

7 Installation and connection

Install the pumps only in accordance with the operating parameters and conditions described in Chapter 4 *Technical data*.

→ Observe the safety instructions (see Chapter 3 *Safety*).



Risk of dangerous gas mixtures during pump operation

Depending on the medium being transferred, breakage of the media-contacting components can result in a dangerous mixture if the medium mixes with the air in the compressor housing or the surroundings.

→ Before using a medium, check the compatibility of the media-contacting components (see 4 *Technical data*) with the medium.

7.1 Installing the pump

→ Store the pump at the same ambient temperature as the installation to mitigate condensation.

Mounting dimensions → For mounting dimensions, see the following figures:

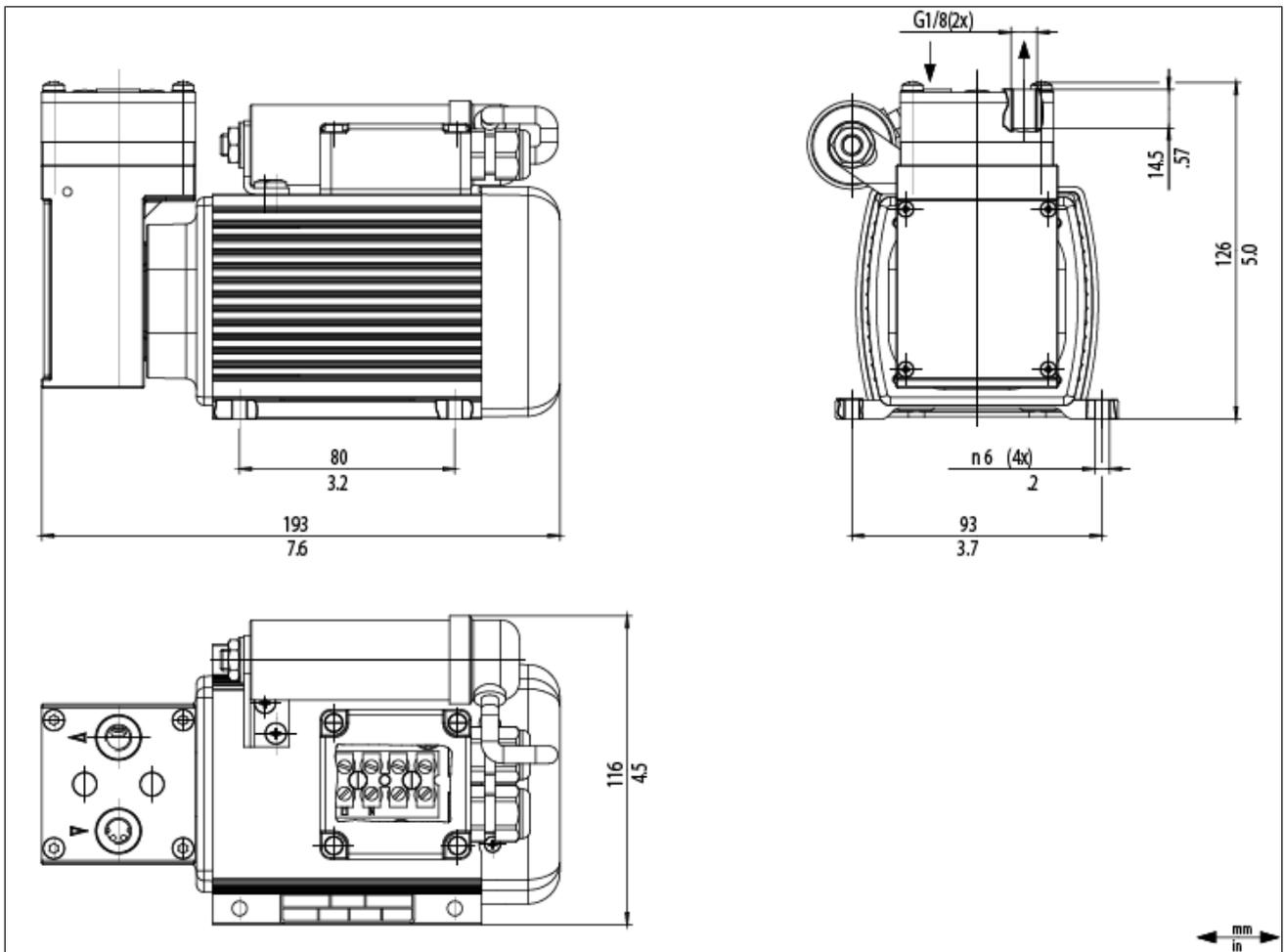


Fig.3 Mounting dimensions N87

Cooling air supply



Danger of burning on hot surfaces

Hot surfaces could occur if the pump overheats.

- When installing the pump, make sure that sufficient cooling air infeed and discharge is ensured.

Immediate environment of the hot pump parts

- During installation, make sure that no combustible or thermally deformable objects are positioned in the immediate area of the hot pump parts (head, drive).

Installation location

- Make sure that the installation location is dry and that the pump is protected from rain, splash water, gushing water, dripping water and other contamination.

- Make sure the pump location in the instrument allows access for periodic maintenance and inspection.



The IP protection class of the pump motor is specified on the type plate.

- Mount the pump at the highest point in the system to prevent condensate from collecting in the pump head.

- Protect the pump from dust.

- Protect the pump from vibration and impact.

7.2 Electrical connection



Danger to life from electric shock

- Only have the pump connected by an authorized specialist.

- Only have the pump connected if the power supply is disconnected.

- When connecting to a power source, observe the applicable regulations, directives, and technical standards.

- Install a device for separating the pump motor from the electrical mains in the electrical installation (e.g., in accordance with EN 60335-1).

- Protect the pump motors, e.g., in accordance with EN 60204-1 (over-current protection, overload protection).



Refer to the type plate for the maximum current consumption of the pump.

- It is recommended that an additional EMERGENCY STOP device be installed.

- Mount the pumps in such a way that it is not possible to touch electrically live parts (electrical connection).

Fastening the connection cables

- Fasten the connection cables so that

- the cables do not come into contact with movable or hot parts.
- the cables cannot be worn or damaged on sharp corners or edges
- no tensile and pressure forces are exerted on the connection point of the cables (strain relief)

Connecting the pump

1. Confirm that the supply voltage meets the parameters listed on the pump nameplate. Refer to the pump nameplate for the rated current consumption.
2. Connect the ground conductor to the pump motor.
3. Connect the motor wires.
4. Open the terminal box cover.
5. For pumps with capacitor motors:
Connect the electrical power cables to terminals L1 and N of the pump motor.
6. Close the terminal box cover again.

7.3 Pneumatic connection



Personal injury or property damage through ejected plugs

If not removed, the plugs on the pressure side of the pump can be ejected during operation by the resulting overpressure.

→ Remove the plugs during installation.

→ Wear appropriate personal protective equipment.

- Connected components → Only connect components to the pump that are designed for the pneumatic data and thermal requirements of the pump. (see Chapter 4 *Technical data*).
- Pump discharge → If the pump is being used as a vacuum pump, safely (relating to the medium and noise) discharge the possibly hot pump discharge via the pneumatic outlet of the pump.

Connecting the pump



Risk of injury from mixing up suction side and pressure side

Mixing up the suction side and pressure side can result in breakage of connected components on the suction side and pressure side.

→ Observe the marking of inlet and outlet on the pump head.

1. Remove the protective plugs from the hose connection threads.
2. Connect the suction line and the pressure line (for mounting dimensions, see Chapter 4 *Technical data*).
3. Lay the suction line and pressure line with a descent so that no condensate can run into the pump.

i Pneumatic noises can be reduced or dissipated by using a silencer (see Chapter 11.2 *Accessories*).

8 Operation

8.1 General



Risk of burns from hot pump parts and/or hot medium

Some pump parts may be hot during or after operation of the pump.

- Allow the pump to cool after operation.
 - Take protective measures to protect against touching hot parts.
-



Injury to eyes

Coming too close to the inlet/outlet of the pump may result in injury to the eyes due to the present vacuum/operating pressure.

- Do not look into the pump inlet/outlet during operation.
-

- Only operate the pumps under the operating parameters and operating conditions as described in Chapter 4 *Technical data*.
- Ensure the proper use of the pumps (See Chapter 2.1 *Proper use*).
- Eliminate the possibility of improper use of the pumps (see Chapter 2.2 *Improper use*).
- Observe the safety instructions (Chapter 3 *Safety*).
- The pumps are built-in devices. Before they are commissioned, it must be ensured that the machines or systems into which the pumps are installed comply with the relevant provisions.



Risk of pump head bursting due to excessive pressure increase

- Do not exceed the maximum permissible operating pressure (see 4 *Technical data*).
- Monitor the pressure during operation.
- If the pressure exceeds the maximum permissible operating pressure of the pump: immediately switch off the pump and remedy the fault (see Chapter 10 *Troubleshooting*).
- Only throttle or regulate the air or gas quantity on the suction line to prevent the maximum permissible operating pressure from being exceeded.
- If the air quantity or gas quantity on the pressure line is throttled or regulated, make sure that the maximum permissible operating pressure at the pump is not exceeded.
- Ensure that the pump outlet is not closed or restricted.

i Excessive pressure, with all of the associated hazards, can be prevented by means of a bypass line with a pressure relief valve between the pressure side and suction side of the pump. Further information is available from KNF Customer Service (contact data: see www.knf.com).



Risk of dangerous gas mixtures during pump operation if diaphragm breaks

If the diaphragm should break, the medium will mix with the air in the compressor housing or in the surroundings.

- Stop pump immediately.
- Replace the diaphragm prior to further operation (see Chapter 9 *Servicing*).

i Operation with open suction-side gas connection can result in contaminants and objects being drawn in.

Pump stoppage → When the pump is at a standstill, establish normal atmospheric pressure in the lines.

Vapors as medium The service life of the diaphragm can be extended, if no condensate forms in the pump. Therefore:

- Perform any work with saturated or near-saturated vapors only with a warm pump.
- KNF recommends: When pumping corrosive media, flush the pump before switching off (see Chapter 9.2.1 *Flushing the pump*) to extend the service life of the diaphragm.

8.2 Information on switching the pump on and off

Switching on the pump

- Ensure that normal atmospheric pressure is present in the lines when switching on.

Switching off/deactivating the pump

- Establish normal atmospheric pressure in the lines (relieve pump pneumatically).
- Recommissioning → Before recommissioning, observe the applicable standards, guidelines, regulations and technical standards at the electrical connection.
- Inspecting the pump → Inspect the pump periodically for external damage or leakage.

9 Servicing



NOTICE

Servicing the pump

Damage to the pumps can result from failure to observe the applicable legal regulations and procedures for the location or intervention by untrained or uninstructed personnel.

- Servicing may only be performed according to the legal regulations (e.g. work safety, environmental protection) and provisions.
- Servicing may only be performed by specialized personnel or trained and instructed personnel.

9.1 Servicing schedule



WARNING

Risk of injury when not using original parts

Failure to use original parts will result in a loss of pump functionality and safety.

The validity of the CE conformity is rendered void if genuine parts are not used.

- Use only genuine spare parts from KNF when performing maintenance/repair work.

Component	Servicing interval
Pump	<ul style="list-style-type: none"> → Inspect the pump periodically for external damage or leakage → Check regularly for conspicuous changes in the noise and vibrations.
Gas connections	<ul style="list-style-type: none"> → Inspect the pump periodically for external damage or leakage

Tab.10

9.2 Cleaning

9.2.1 Flushing the pump

When transferring dangerous and environmentally hazardous media, KNF recommends flushing the pump with air at atmospheric pressure for a few minutes prior to switch-off (if necessary for safety reasons: with an inert gas) to extend the service life of the diaphragm.

- Discharge the media safely.

9.2.2 Cleaning the pump



CAUTION

Risk of burns from hot pump parts

The pump head or motor may still be hot after operation of the pump.

→ Allow the pump to cool after operation.



WARNING

Health hazard due to dangerous substances in the pump

Depending on the medium being transferred, caustic burns or poisoning is possible.

→ Wear protective equipment if necessary, e.g., protective gloves, goggles.

→ Clean the pump with suitable measures.



NOTICE

During cleaning work, ensure that no fluids enter the interior of the housing.

→ Only clean the pump with a dry wiping cloth. When cleaning, use no solvents if possible as these can affect the plastic parts.

→ Only use solvents during cleaning if head materials are not affected (ensure resistance of the material).

→ If compressed air is present, blow out the components.

9.3 Replacing the structured diaphragm and valve plates/ seals (TT design)

Requirements → Disconnect the motor from mains and ensure that it is voltage-free.

→ Clean the pump and free the pump of hazardous materials.

Spare parts

Spare part*	Item designation**	Number per pump head
Structured diaphragm	(7) in Fig. 4	1
Valve plates/seals	(12) in Fig. 4	2

Tab.11 *According to spare parts list, Chapter 11.1 Spare parts

Tool and material

Quantity	Tool/material
1	TORX® screwdriver T10
1	TORX® screwdriver T20
1	Phillips screwdriver
	Thread locker (Loctite 222 or comparable product)
1	Felt-tip pen

Tab.12

Information on the procedure → Always replace diaphragm and valve plate/seals together to maintain the performance of the pump.




WARNING

Health hazard due to dangerous substances in the pump

Depending on the medium being transferred, caustic burns or poisoning is possible.

→ Wear protective equipment if necessary, e.g., protective gloves, goggles.


→ Clean the pump with suitable measures.

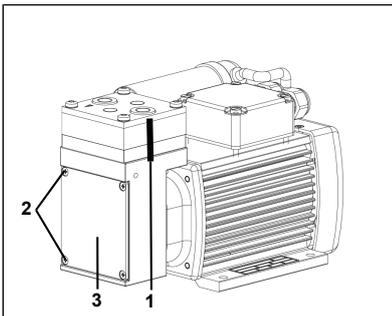

Removing pump head

1. Mark the pressure plate, head plate, intermediate plate and housing with a felt-tip pen (Fig. /1) to ensure proper mounting.
2. Remove the housing cover (Fig. /3).
3. Loosen the four fastening screws (Fig. 4/2) and remove the housing cover (Fig. 4/3).

The counterweight is visible.

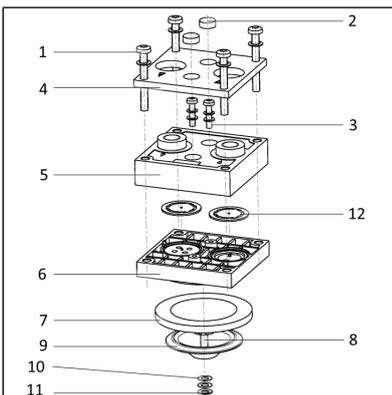

4. Loosen the four screws (Fig. 4/1) and remove the pressure plate (Fig. 4/4).
5. Remove the screw covers (Fig. 4/2); loosen the two screws (Fig. 4/3) and remove the head plate (Fig. 4/5) and the intermediate plate (Fig. 4/6).

The structured diaphragm (7) is visible.

Replacing the structured diaphragm

1. Turn the pump to the side; this keeps the shim rings (10) and the disc springs (11) from falling into the pump chamber when removing the structured diaphragm (7).
2. Move the structured diaphragm (7) by turning the counterweight to the upper reversal point.
3. Lift the structured diaphragm (7) onto opposing side edges; hold the structured diaphragm and unscrew it counterclockwise.
4. Remove the diaphragm support (9), shim ring(s) (10) and disc spring (11) from the threaded bolt (8) of the structured diaphragm and keep in a safe place.
5. Check all parts for soiling and clean them if necessary.

Fig.4 Replacing the structured diaphragm

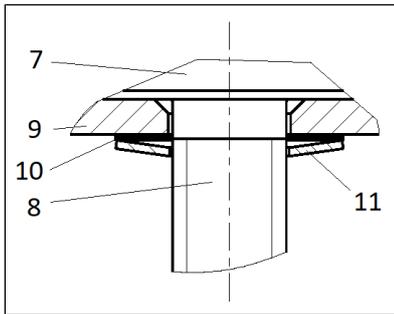


Fig.5 Aligning disc spring

6. Push diaphragm support (9), shim ring(s) (10) and disc spring (11) onto the threaded bolt (8) of the new structured diaphragm (7) in this order (see Fig. 5).

i The edge of the disc of the disc spring (11) must be aligned with the structured diaphragm (7).

7. Move the connecting rod (connection part between drive shaft and structured diaphragm) to the upper reversal point by turning the counterweight.
8. Screw the structured diaphragm with diaphragm support, shim ring(s) and disc spring clockwise onto the connecting rod and hand tighten.

Replacing valve plates/seals

1. Remove the valve plates/seals (12) from the intermediate plate (6).
2. Check the valve seats, intermediate plate (6) and head plate (5) for soiling and damage. Clean the parts if necessary.
3. Contact KNF in the event of unevenness, scratches or corrosion. Order and replace damaged parts.
4. Insert the new valve plates/seals (12) in the valve seats of the intermediate plates (6).
5. Place the head plate (5) on the intermediate plate according to the felt-tip pen marking.
6. Check the centering of the head plate (5) by means of a slight lateral movement.
7. Connect head plate (5) and intermediate plate (6). To do this, screw in screws (3; Fig. 6: A + B) with disc spring (11) and shim rings (10) until the screw head is seated flat, i.e., until the first resistance is encountered. Apply a small amount of thread locker (Loctite 222 or comparable product) to the threads of the screws.

Mounting the pump head

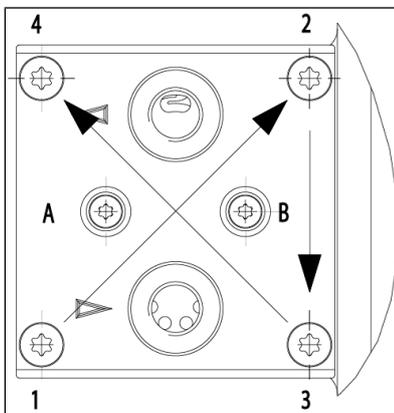


Fig.6 Tightening cap head screws

1. Place the pump head on the housing according to the felt-tip pen marking.
2. Place the pressure plate (4) on the head plate according to the felt-tip pen marking.
3. Screw in the screws (1) with disc spring (11) and slightly tighten them crosswise in the order 1-2-3-4 (see Fig. 6) with 1 Nm.
4. Check for ease of movement of the pump by turning the counterweight.
5. Tighten the screws (1) with disc spring (11) crosswise in the order 1-2-3-4 (see Fig. 6). Tightening torque: 2.1 Nm.
6. Then tighten the screws (3; Fig. 6: A + B) another quarter turn (tightening torque: 30 Ncm) and mount the screw covers (2).

Final steps

1. Mount the housing cover.
2. Check the pump head and pneumatic connections for leaks.
3. Properly dispose of the replaced structured diaphragm and valve plates/seals.

i To ensure the required gas tightness of the pump following servicing, a leak test is to be performed.



Risk of injury and poisoning from leaks

→ Before recommissioning the pump, check the pump heads and pneumatic connections for leaks. Leaks may lead to poisoning, chemical burns or similar injuries.

9.4 Changing the structured diaphragm and valve plates/ seals (S_-/A_ design based on project)

- Requirements → Disconnect the motor from mains and ensure that it is voltage-free.
 → Clean the pump and free the pump of hazardous materials.

Spare parts

Spare part*	Item designation**	Number per pump head
Structured diaphragm	(7) in Fig. 7	1
Valve plates/seals	(12) in Fig. 7	2

Tab.13 *According to spare parts list, Chapter 11.1 Spare parts

Tool and material

Quantity	Tool/material
1	TORX® screwdriver T20
1	Phillips screwdriver
	Thread locker (Loctite 222 or comparable product)
1	Felt-tip pen

Tab.14

- Information on the procedure → Always replace diaphragm and valve plate/seals together to maintain the performance of the pump.



Health hazard due to dangerous substances in the pump

Depending on the medium being transferred, caustic burns or poisoning is possible.

- Wear protective equipment if necessary, e.g., protective gloves, goggles.
- Clean the pump with suitable measures.

Removing the pump head

1. Mark the head plate, intermediate plate and housing with a felt-tip pen (*Changing the structured diaphragm and valve plates/seals (S_-/A_ design based on project)*/1) to ensure proper mounting.

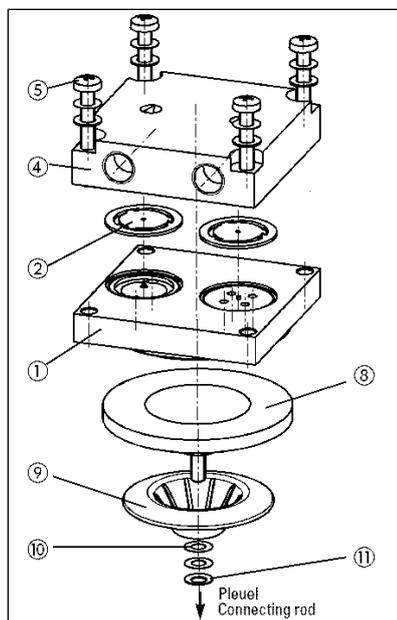


Fig.7 Replacing the structured diaphragm

2. Remove the housing cover (*Changing the structured diaphragm and valve plates/seals (S_-/A_ design based on project)*3).
3. Loosen the four fastening screws (Fig. 4/2) and remove the housing cover (Fig. 4/3).

The counterweight is visible.

4. Loosen the four screws (Fig. 7/5).
5. Remove the head plate (Fig. 7/4) and the intermediate plate (Fig. 7/1).

The structured diaphragm (8) is visible.

Replacing the structured diaphragm

1. Turn the pump to the side; this keeps the shim rings (10) and the disk springs (11) from falling into the pump chamber when removing the structured diaphragm (8).
2. Move the structured diaphragm (8) by turning the counterweight to top dead center.
3. Lift the structured diaphragm (8) onto opposing side edges; hold the structured diaphragm and unscrew it counterclockwise.
4. Remove the diaphragm support (9), shim ring(s) (10) and disk spring (11) from the threaded bolt of the structured diaphragm (8) and keep in a safe place.
5. Check all parts for contamination and clean them if necessary.
6. Push the diaphragm support (9), shim ring(s) (10) and disk spring (11) onto the threaded bolt of the new structured diaphragm (8) in this order (see *Changing the structured diaphragm and valve plates/seals (S_-/A_ design based on project)*).

i The edge of the disc of the disk spring (11) must be aligned with the structured diaphragm (8).

7. Move the connecting rod (connection part between drive shaft and structured diaphragm) to top dead center by turning the counterweight.
8. Screw the structured diaphragm with diaphragm support, shim ring(s) and disk spring clockwise onto the connecting rod and hand tighten.

Replacing the valve plates/seals

1. Separate the head plate (4) from the intermediate plate (1).
2. Remove the valve plates/seals (2) from the intermediate plate (1).
3. Check the valve seats, intermediate plate (1) and head plate (4) for contamination and damage. Clean the parts if necessary.
4. Contact KNF in the event of unevenness, scratches or corrosion. Order and replace damaged parts.

i The valve plates/seals are identical for the pressure and suction side. The same applies for the top and bottom of the valve plates/seals.

5. Insert the new valve plates/seals (2) in the valve seats of the intermediate plates (1).
6. Place the head plate (4) on the intermediate plate according to the felt-tip pen marking.
7. Check the centering of the head plate (4) by means of a slight lateral movement.

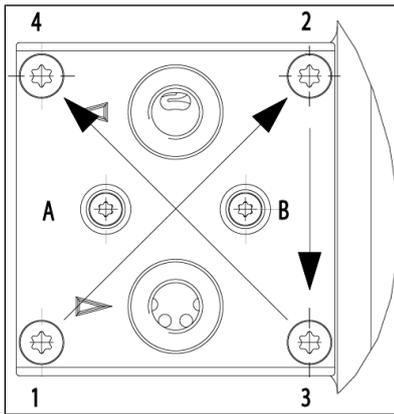


Fig.8 Tighten cap head screws
(head plate shown as an example
(screws (A) and (B) only in N87TTE
design))

Mounting the pump head

1. Place the pump head on the housing according to the felt-tip pen marking.
2. Screw in the screws (5) with disk spring and slightly tighten them crosswise in the order 1-2-3-4 (see Fig. 8) with 1 Nm.
3. Check for ease of movement of the pump by turning the counterweight.
4. Tighten the screws (5) with disk spring crosswise in the order 1-2-3-4 (see Fig. 8). Tightening torque: 2.1 Nm.

Final steps

1. Mount the housing cover.
2. Check the pump head and pneumatic connections for leaks.
3. Properly dispose of the replaced structured diaphragm and valve plates/seals.



To ensure the required gas tightness of the pump following servicing, a leak test is to be performed.



WARNING

Risk of injury and poisoning from leaks

- Before recommissioning the pump, check the pump heads and pneumatic connections for leaks. Leaks may lead to poisoning, chemical burns or similar injuries.

10 Troubleshooting



Danger: electric shock can be life-threatening.

- All work on the pump may only be performed by an authorized specialist.
- Before working on the pump: Disconnect the pump from the power supply.
- Check and ensure that no voltage is present.

→ Allow the pump to cool before troubleshooting.

→ Check the pump (see following tables).

Unit not pumping	
Cause	Troubleshooting
Pump is not connected to the electrical mains.	→ Connect the pump to the electrical mains.
No voltage in the electrical mains.	→ Check the circuit breaker for the room and switch it on if necessary.
Pneumatic connections or lines are blocked.	→ Check the connections and lines. → Remove the blockage.
External valve is closed or filter is plugged.	→ Check external valves and filters.
Condensate has collected in the pump head.	→ Separate the source of the condensate from the pump. → Flush the pump with air at atmospheric pressure for a few minutes (if necessary for safety reasons: with an inert gas). → Install the pump at the highest location in the system.
Diaphragm or reed valves are worn or defective.	→ Change the diaphragm and valve springs (see Chapter Changing diaphragm and reed valves).

Tab.15

Flow rate, pressure or vacuum too low	
The pump does not reach the performance stated in the technical data or data sheet.	
Cause	Troubleshooting
Condensate has collected in the pump head.	<ul style="list-style-type: none"> → Separate the source of the condensate from the pump. → Flush the pump with air at atmospheric pressure for a few minutes (if necessary for safety reasons: with an inert gas). → Install the pump at the highest location in the system.
There is overpressure on the pressure side and at the same time vacuum or pressure above atmospheric pressure on the suction side.	<ul style="list-style-type: none"> → Change the pneumatic conditions.
Pneumatic lines or connection parts have insufficient cross section or are throttled.	<ul style="list-style-type: none"> → Disconnect the pump from the system to determine the output values. → Eliminate any throttling (e.g., valve) if necessary. → Use lines or connection parts with a larger cross section if necessary.
Leaks occur at pneumatic connections, lines or pump head.	<ul style="list-style-type: none"> → Eliminate the leaks.
Pneumatic connections or lines are completely or partially clogged.	<ul style="list-style-type: none"> → Check the pneumatic connections and lines. → Remove any parts or particles that are causing blockages.
Head parts are soiled.	<ul style="list-style-type: none"> → Clean the head components.
Operating diaphragm broken	<ul style="list-style-type: none"> → Shut down the pump immediately.
Diaphragm or reed valves are worn or defective.	<ul style="list-style-type: none"> → Change the diaphragm and valve springs (see Chapter Changing diaphragm and reed valves).

Tab.16

Pump exhibiting changed running noises and vibrations.	
Cause	Troubleshooting
Pump bearing worn or defective.	<ul style="list-style-type: none"> → Determine the cause. → Contact KNF Customer Service.
Drive worn or defective.	<ul style="list-style-type: none"> → See operating instructions for the motor.

Tab.17

Fault cannot be rectified

If you are unable to identify any of the specified causes, send the pump to KNF Customer Service (contact data: see www.knf.com).

1. Flush the pump with air for a few minutes (if necessary for safety reasons: with inert gas) at atmospheric pressure to free the pump head of dangerous or aggressive gases (see Chapter 9.2.1 *Flushing the pump*).
2. Clean the pump (see Chapter 9.2.2 *Cleaning the pump*).
3. Send the pump together with completed Health and Safety Clearance and Decontamination Form to KNF, stating the nature of the transferred medium.

11 Spare parts and accessories

i To order spare parts and accessories, please contact your KNF sales partner or KNF Customer Service (contact data: see www.knf.com).

11.1 Spare parts

Spare part	Order number
Valve plate/seal	057172
Structured diaphragm	044973

Tab.18 N87TTE spare parts

11.2 Accessories

Description	Order number
Hose connector for hose ID 6 (PVDF, G 1/8)	014052
Male connection OD 6 (PVDF, G 1/8)	014049

Tab.19

12 Returns

Preparing for return

1. Flush the pump with air for a few minutes (if necessary for safety reasons: with inert gas) at atmospheric pressure to free the pump head of dangerous or aggressive gases (see Chapter 9.2.1 *Flushing the pump*).

i Please contact your KNF sales partner if the pump cannot be flushed due to damage.

2. Remove the pump.
3. Clean the pump (see Chapter 9.2.2 *Cleaning the pump*).
4. Send the pump together with the completed Health and Safety Clearance and Decontamination Form to KNF, stating the nature of the transferred medium.
5. Pack the device securely to prevent further damage to the product. If necessary, request original packaging for a fee.

Returns

KNF shall undertake to repair the pump only under the condition that the customer presents a certificate regarding the medium that is pumped and the cleaning of the pump. Please follow the instructions at knf.com/repairs here.

Contact your KNF sales partner directly if you require additional support for your return service.

KNF worldwide

You can find our local KNF partners at: www.knf.com